

# **200 eV to 30 keV** Medium Energy, Uniform Flood Electron Beam

# FOR USE IN:

- Spacecraft Charge-up Studies
- Surface Charging Studies
- Beta Decay Stimulation
- Surface Physics Studies
- Vacuum Physics Experiments

# FEATURES / OPTIONS:

- Energy Range: 200 eV to 30 keV
- Beam Currents to 1 mA
- ➢ Wide Angle Uniform Beam
- Divergence Control
- Rotatable 2.75" CF Mounting
- User-Replaceable Firing Units
- Computer / Remote Control
- Custom Apertures



EGF-4104 Electron Flood Gun (shown with deflection option)

The Kimball Physics EGF-4104 Electron Gun, with its matching EGPS-4104 Power Supply, is intended for use in a variety of UHV charging, space physics, vacuum physics, surface physics, and nuclear simulation applications. It is a complete subsystem ready to attach and turn on. Maximum flexibility is achieved in a minimum of space; the entire unit mounts through a single standard 2.75" inch CF port.

Electrons are generated at negative high potential, and the user's target is typically set at ground potential. Both beam energy and beam current are adjustable over wide ranges. The gun uses a space-charge-limited refractorymetal cathode to generate a uniform flood beam and the design allows generation of the beam down to low energies, and very low currents. A high current option provides beam currents up to 1 mA. Beam divergence is partially controllable electronically over the full range of the electron energy. Optional cathodes include barium oxide discs (BaO, low light, low energy spread) or yttriacoated iridium discs (Y2O3 - Ir, rugged, may survive brief loss of vacuum). Except for barium oxide, cathodes are not damaged by repeated exposure to atmospheric gases or water vapor when cold. Cathode lifetime is a function of vacuum conditions and beam current as related to cathode temperature. Cathode lifetime at low currents in good vacuum may be in the many hundreds of hours, or even over a thousand hours.

UHV technology is used throughout. The gun may be completely disassembled for cleaning and repair. The cathode firing unit assembly (which includes the cathode, cathode mount, apertures, and Wehnelt) is user-replaceable; and the firing unit assemblies may be sent back to Kimball Physics for rebuilding. The gun can be run in vacuums from10<sup>-11</sup> torr to 10<sup>-5</sup> torr. The gun may be baked up to 350°C with cables removed (150°C with magnetic deflection option). Non-standard mountings are available. Various designs of stand-alone Faraday cups are available.

The EGPS-4104 Power Supply contains all power supplies necessary to generate the required voltages to run an EGF-4104 Electron Flood Gun, including the Energy, Focus, Source, Grid, Anode, and optional Deflection supplies. All power supplies are electronically regulated. A computer control option allows for control of all potentials from an interface at ground potential, via 0 to 10 V analog inputs.

The FlexPanel provides a digital display screen and a keypad controller for programming control on the front panel. Rear panel connectors allow remote /computer control and metering of all gun power supplies. A mini-USB and RS-232 serial ports, and an analog input/output connector are included on standard power supply units. All common computer interface bus types can be accommodated, by use of appropriate digital to analog converters. RS-422/485 conversion is possible.

An optional LabVIEW<sup>TM</sup> computer program designed for the EGF-4104 is available for remote computer control and metering. Software is available in two options: 1) using National Instrument DAQ modules and the 50-pin connector on the EGPS-4104, or 2) via a mini-USB or simple serial connector interface. The program provides a virtual panel of controls and real-time metering on the user's computer screen.



EGPS-4104B Electron Gun Power Supply with FlexPanel controller.

| EGF-4104 ELECTRON GUN SPECIFICATIONS  |   |
|---|---|
| BEAM ENERGY   | 200 eV to 30 keV (Independently adjustable)   |
| BEAM CURRENT  | Standard: 1 nA to 100 µA<br>(Independently adjustable) High<br>current option: 10 nA to 1 mA*   |
| ENERGY SPREAD   | Approx. cathode thermal spread,<br>calculated<br>Ta - 0.5eV<br>Y2O3 - 0.4eV BaO - 0.3eV   |
| SPOT SIZE   | Typical - 15 to 450+ mm   |
| WORKING DISTANCE  | Typical - 100 to 1000 mm  |
| BEAM DEFLECTION   | Optional: Magnetic quadrupole<br>(outside vacuum)   |
| PULSE CAPABILITY<br>(using appropriate<br>pulse generator,<br>not included) | Dual Grid Power Supply: pulse width ~2 µs to DC rise/ fall ~500 ns, rep rate to 5 kHz (TTL required)  |
| BEAM UNIFORMITY   | Variable. Optional beam washing or mask aperture  |
| FIRING UNIT   | User-replaceable Firing Unit Cartridge<br>includes cathode and Wehnelt/Grid (G-<br>1) assembly  |
| CATHODE TYPE  | Standard: Tantalum<br>Optional: Low-light Barium Oxide<br>Yttria-coated Iridium   |
| BEAM<br>ALIGNMENT   | Mechanical alignment with internal firing unit alignment. Optional mechanical alignment with $a \pm 2^{\circ}$ Port Aligner.  |
| MOUNTING  | 2 <sup>3</sup> / <sub>4</sub> inch rotatable CF, including both tapped and clear mounting holes   |
| INSERTION LENGTH  | 0 mm  |
| GUN DIMENSIONS  | 85 mm OD (excluding Lemo<br>connector) x 205 mm   |
| FEEDTHROUGHS  | Multi-pin brazed ceramic, threaded stainless steel shell  |
| CABLES /<br>CONNECTORS  | Multi-conductor high voltage fully<br>ground-shielded cable with mating<br>aluminum shell connector, to connect<br>gun and power supply. Standard length:<br>3 m, Optional: 5 m |
| MAXIMUM BAKEOUT   | 350°C with cables removed<br>150oC with magnetic deflection option  |

| EGPS-4104 POWER SUPPLY SPECIFICATIONS  |  |
|--|--|
| OUTPUT                                 | All necessary voltages to drive the<br>EGF-4104 Electron Gun (in<br>combination with H.V. Power<br>Supply*)  |
| ENERGY SUPPLY<br>STABILITY             | $\pm 0.01\%$ per hour $\pm 0.02\%$ per 8 hours at full output  |
| BEAM STABILITY                         | $\pm 0.1\%$ per hour with Emission Current<br>Control (ECC) or $\pm 10\%$ per hour<br>without ECC  |
| CONTROLS                               | FlexPanel controls: Energy, Source,<br>Grid (G-1), 1st Anode, Focus,<br>Emission Current Control (ECC),<br>Optional X and Y Deflection.  |
| METERING                               | FlexPanel digital meters: Energy,<br>Source Voltage, Source Current, Grid<br>Voltage, Anode Voltage, Focus<br>Voltage, Emission Current, Optional X<br>and Y Deflection          |
| COMPUTER/REMOTE<br>CONTROL & METER     | Controls: 0 to +10V (-10V to +10V,<br>for deflection) Metering: 0 to +2V (-<br>2V to +2V, for deflection) Toggle<br>switches: 0 or +5V   |
| SOFTWARE                               | Standard configuration designed for<br>serial USB or RS-232 connections.<br>Optional: National Instruments<br>LabVIEW <sup>TM</sup> file, designed to run with<br>NI DAQ modules |
| INPUT                                  | 115 VAC or 230 VAC, 50 to 60 Hz<br>single phase, 250 VA  |
| ENVIRONMENT                            | Temperature: 0 to 40°C, Relative<br>humidity: 0 to 75% RH non-<br>condensing, Classified as a pollution<br>degree 2, installation category                                       |
| DIMENSIONS<br>(width x height x depth) | EGPS-4104: 17 in. x 7 in. x 22 in.<br>excluding handles (432 mm x 178 mm<br>x 560 mm); 19 in.(495 mm) rack<br>mountable.   |
| WEIGHT                                 | Appx. 35 lbs. (16 kg)  |

### OPTIONAL HARDWARE RASTER SPECIFICATIONS

| RASTER<br>GENERATOR | Frequency in X and Y directions can be<br>independently set. X freq. is up to 500 Hz, Y<br>freq. 100Hz. When X freq. exceeds 100Hz the<br>raster angle is reduced by 25%.<br>All parameters controllable via serial USB or<br>RS-232, RS-485, analog input, or computer<br>control with LabVIEWTM software option. |
|---------------------|--|
|                     |  |



4



EGF-4104 ELECTRON GUN with optional deflection

# References

For more information on electron sources / gun operations (and the technical bulletins and additional documents listed below), please visit the Resources and Documents section of our website.

**General Operating Hints** 

Operating Instructions, Typical LabVIEW<sup>™</sup> Electron Gun Systems

**Beam Pulsing Options** 

Note: A comprehensive custom manual is supplied with each system.

#### Notes:

5

- 1. Charts /graphs show typical performance, data is for guidance only
- 2. It is not necessarily possible to achieve all maximum specifications simultaneously.
- 3. Specifications Subject to Change Without Notice.
- 4. DE Altobelli, DT Taylor 04/15/2025

Document EGF-4104\_EGPS-4104\_2025\_0415 COPYRIGHT KIMBALL PHYSICS 2025, ALL RIGHTS RESERVED