


**KIMBALL PHYSICS**   
*Excellence in Electron and Ion Optics*

**KIMBALL PHYSICS ELECTRON GUNS**  
TUTORIAL AND RESOURCE

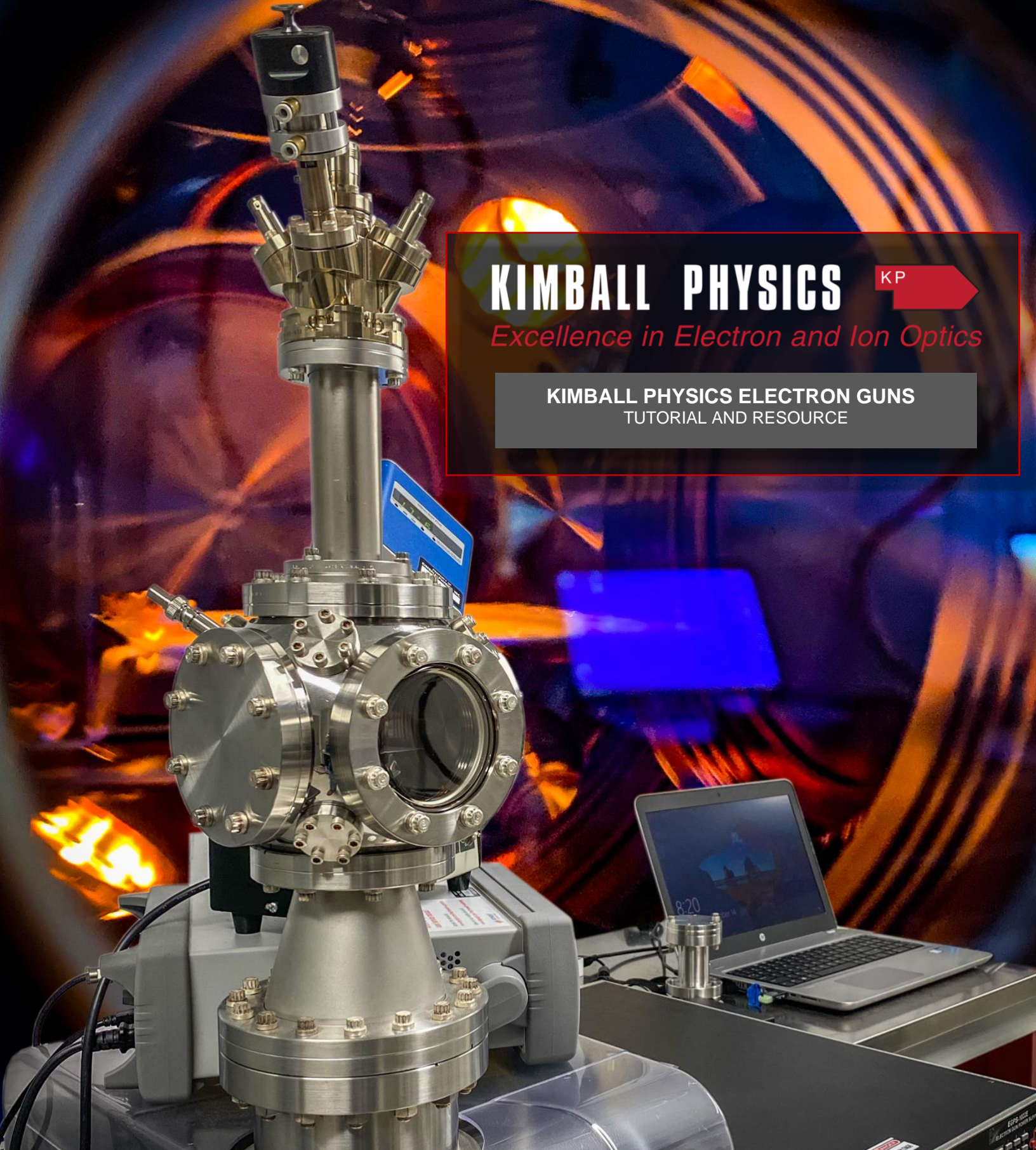

---

## Table of Contents

Introduction	3
Basic Components Electron Gun System	3
Electron Guns: Basic Anatomy	4
Triode	
Electron Emitter	
Electrodes and Electric Fields	
Anode and Cathode	
Wehnelt Cylinder or Grid	
Electron Optics Overview	7
Supporting Systems	7
Physical Systems	
Vacuum Systems Overview	
Power Systems Overview	
Detectors	
Electron and Ion Gun Systems (201)	10
Beam Energy	
Emission and Beam Current	
Beam Spot Size	
Working Distance	
Beam Divergence	
Beam Uniformity and Current Density	
Brightness	
Electron Optics	12
Beam Focusing	
Zoom Lens	
Deflection /Rastering	
Magnetostatic Lens	
Cross Over Points	
Beam Operation and Performance Factors	15
Beam On / Off	
Grid Cut Off	
Beam Pulsing	
Electron Sources	17
Other Types of Electron Emission	
Photoelectron Emission	
Field Emission	

Ion Sources and Systems	18
Gas Ionization	
Plasma	
Solid Ionization	
Ion Beam	
Frequently Asked Questions:	
Gun Operations	20
How is the Beam Moved on the Target?	
Mechanical Alignment	
Deflection	
Rastering	
Beam Rocking	
How is the Electron or Ion Beam Observed?	
Phosphor Screen	
Faraday Cup	
How are the Power Supplies Controlled?	
Manual or Remote Control	
Feedback Control (ECC)	
How is the Gun Connected to a Vacuum System?	
Mounted Guns	
Unmounted Guns	
How are Guns Tested?	
Data	
References	24
Glossary	26
Notes	36

---

## Electron Gun Tutorial and Resource

### Introduction

Kimball Physics specializes in the design and manufacturing of precision high-tech scientific instruments with over 50 years of experience in ultra-high-vacuum electron and ion optics. Our expertise is in high stability electron emitters, precision electrostatic and magnetostatic optics, along with state-of-the-art vacuum chambers and fittings. Electron and ion systems are optimized for beam energies ranging from 1 eV to 100 keV and beam currents from femtoamperes to amperes.

Our goal with this informal resource is to provide a fundamental knowledge base for understanding the basic terminology, principles and parameters, and specifications of electron and ion beam systems.

Many scientists, engineers and graduate students would like to incorporate electron and ion beam systems into applications within their specific fields and need to “get up to speed” quickly with this fundamental information. With this information resource, our intent is to help you move more quickly towards your solution and to choose the electron or ion system and components that would work best for your application.

The first section, Introduction to Electron Beam Systems (101), is a ground level introduction to the terminology and basic mechanisms related

to electron beam systems. The next section, Electron and Ion Gun Systems (201), provides more advanced details about the electron and ion beam parameters and the associated instrumentation. The intent of this section is to provide more specific and detailed information about the systems to enable the interpretation of the technical specification documents available for our electron and ion gun products and options. The last section includes basic Gun Operations and Questions. Advanced Topics (301) will be added over time and will dive into more detail about the scientific and technical aspects of chosen topics related to the systems.

Please excuse any errors or areas that are not complete; this document is always a work in progress!

### Basic Components of an Electron Gun System

Let's get started with a dissection of an electron gun system with basic description of how the various elements work and come together. For this first pass through the information, the description of each element will be basic. Later in the document, we will add more details, parameters, and technical description.

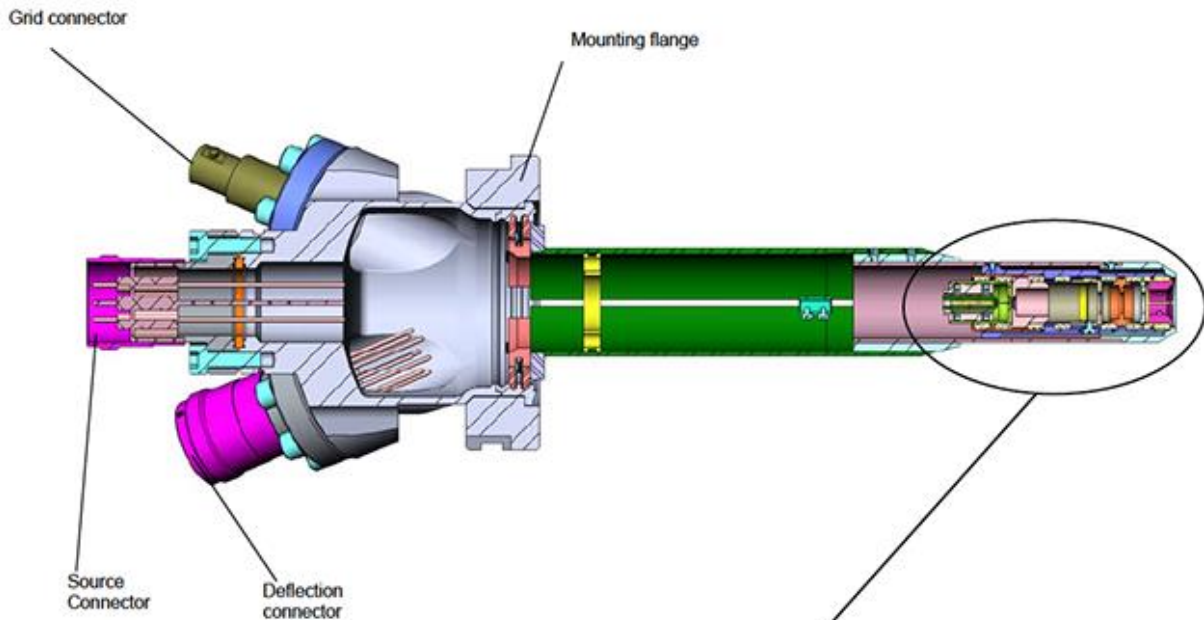


Figure 1. Example of a Kimball Physics Electron Gun.


## Electron Guns: Basic Anatomy

The figures included will provide an overview of a typical electron gun. The first figure above (Figure 1), is an exhibit of a typical electron gun. Figure 2 below is a 3D model, with a cross-section of a typical electron gun with the

elements labeled. Figure 3 below is a simplified schematic of the electron gun components and applied voltages. Figure 4 shows a schematic block diagram of the typical power supplies (represented as circles) and how the voltages are typically applied to the parts of the gun to produce and control the electron beam. Please



**Figure 2.** 3D solid model with a zoomed in area with more detailed three-dimensional drawing showing a labeled cross-section of a typical electron gun components.

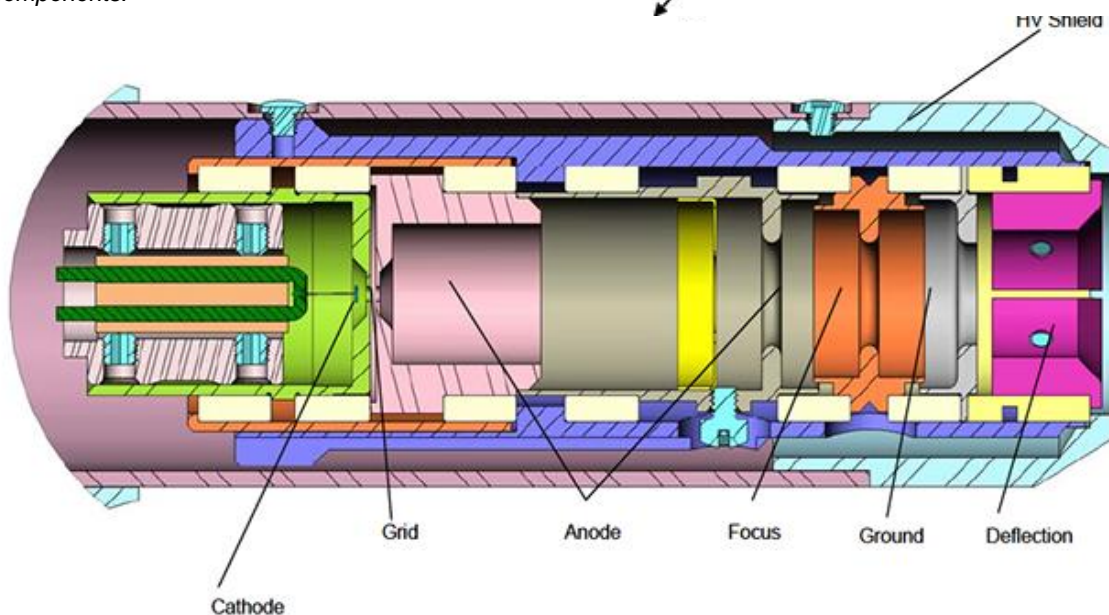

reference these figures as we progress through the discussions that follow.

### Triode

The electron gun is a descendant of the early vacuum tube, and some parts still utilize the same names although they are somewhat different in structure.

In a triode vacuum tube, which is the basis for many electron guns, there are three electrodes- 1) the *cathode* that generates free electrons, 2) the *anode* which attracts and accelerates the electrons to a displaced point in the beam direction, and 3) a *grid* (in electron guns know as a Wehnelt grid) to provide control of the electron emission from the cathode.

### Electron Emitter

The simplest source of the electrons in an electron gun cathode could be a metallic wire or filament such as tungsten bent into a hairpin shape, where we apply a small voltage to pass a current through the wire's resistance to heat the wire.

Depending on the temperature and characteristics of the wire, such as its work


function (we talk more about work function, cathode materials, designs, and parameters in our cathode tutorial document), some of the electrons in the wire will have enough energy to escape their outer valance shells, become free electrons and then possibly emerge to the surface of the cathode and into the partial vacuum space just surrounding it (*thermionic emission*). The electrons initially exit from the material surface in several random directions and at low energies.

### Electrodes and Electric Fields

Electrodes are simply conductors with geometries that interface with non-conductors in the system (such as the partial vacuum space). They are typically used to establish a point of voltage potential and in this case, to create electric field gradients within the electron gun's physical configuration. In addition to the cathode, the anode and Wehnelt grid, there are typically several additional electrodes present in the system that modify the path of the electron beam.

### Anode and Cathode

To facilitate the extraction of electrons emitted from the cathode, and then organize and

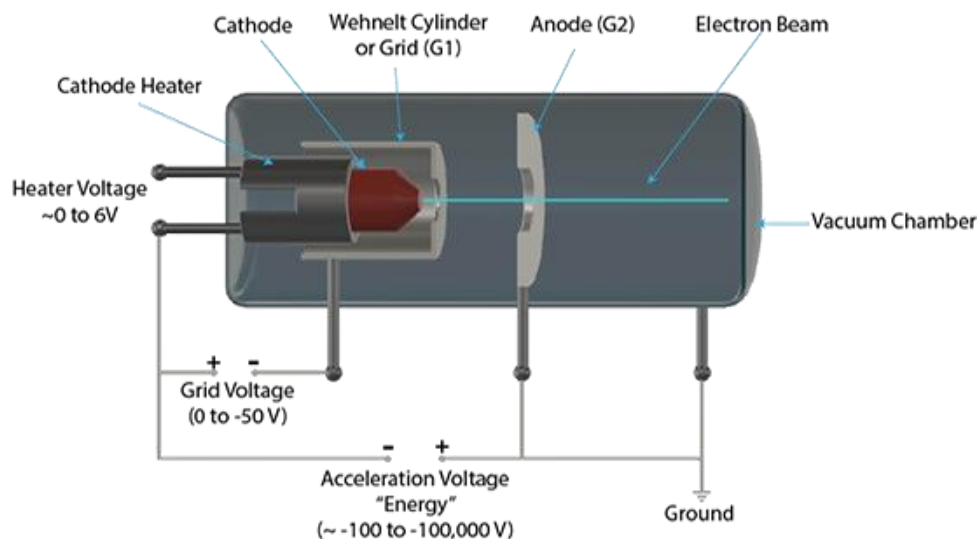

**Figure 3.** Schematic view of simplified electron gun with labelled components.


accelerate them into a beam with more uniform characteristics, a strong primary electric field is created in the region.

The electron source or cathode (the negative electrode) is typically held at a large negative potential (such as -1000 V or more) with respect to the anode (the positive electrode), where the anode is in a position at some distance and in the intended beam direction. The anode could be in the shape of a conductive disk (if it was to be the target) or ring geometry if the beam was to pass through or by it.

The anode's potential is often held at ground (0 V), which is at a much higher potential than the cathode's negative potential. As the electrons are emitted from the source and accelerated

towards the anode by the electric field established between the cathode and anode, they become somewhat more organized into a beam like structure. As the beam continues past the anode, other electron optics techniques and mechanisms are implemented to focus and modify the beam trajectory and other beam characteristics. The potential difference between the cathode and anode largely determines the overall energy of the electron beam.

The Schottky effect or field enhanced thermionic emission is a phenomenon related to how the electric field magnitude lowers the surface barrier for electron emission and allows increased emission current at a given temperature.

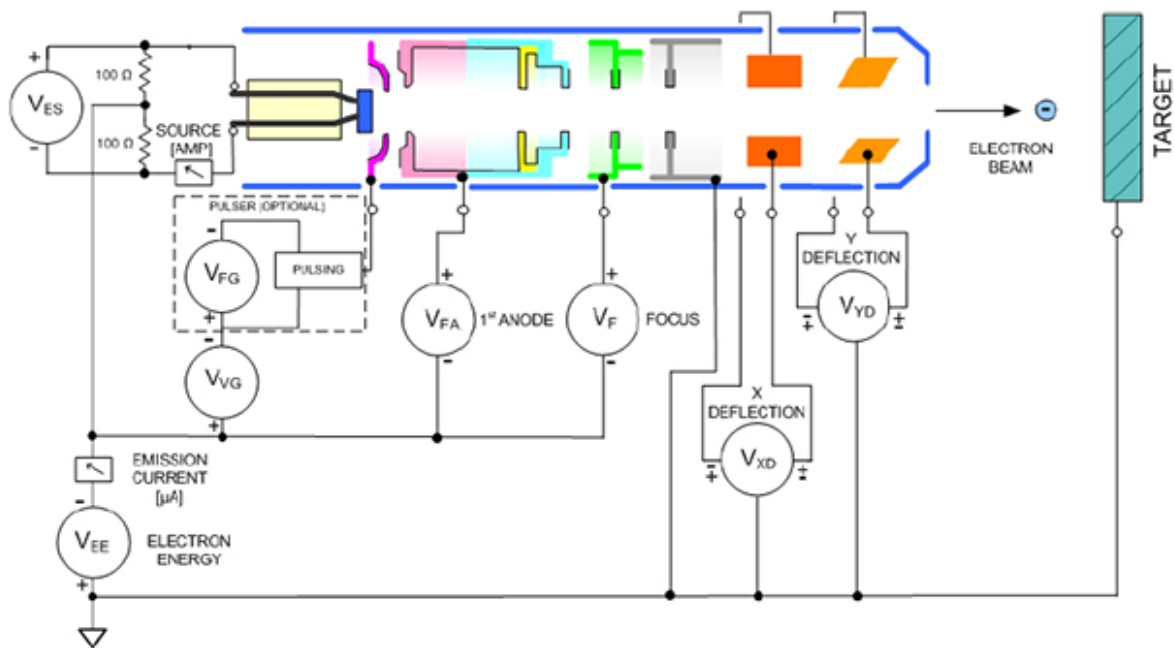

Voltage Ranges	
V <sub>EE</sub> ENERGY (HIGH)	up to -2 keV
(LO)	up to -200 eV
V <sub>VG</sub> GRID	0 to -50 V
V <sub>FA</sub> 1 <sup>ST</sup> ANODE	0 to 200 V
V <sub>F</sub> FOCUS	0 to 2000 V
V <sub>ES</sub> SOURCE	0 to 3V @ 3A
V <sub>XD</sub> DEFLECTION	-150 V to +150 V
V <sub>YD</sub> DEFLECTION	-150 V to +150 V
FIXED GRID (OPTIONAL)	-50V

Figure 4. Simplified schematic of electron gun with typical relative potentials between components.

---

### Wehnelt Cylinder or Grid

To help focus and control the release of electrons from the emitting source, a cup-shaped-cylindrical element with a hole or aperture in its center is placed in close proximity to the cathode (emitting) surface. This is commonly referred to as the Wehnelt cylinder or grid, and it is typically kept at a potential that is even more negative than the adjacent cathode source.

The potential difference between the cathode electron source and the Wehnelt cylinder creates another, more localized electric field, which is somewhat opposite in direction when compared to the primary electric field between the cathode electron source and the more distant anode.

At low magnitude, this electric field will suppress emission from the lateral surfaces of the electron source to initially help focus the random electrons into a more beam-like configuration. The electric field (formed between the cathode and Wehnelt) can also be increased to fully suppress emission of electrons from the electron source and can be used to create a pulsed electron beam. The Wehnelt or grid is the first element in a typical series of devices to control the electron beam that has been formed.

### Electron Optics Overview

An interesting property of a glass lens is its ability to bend electromagnetic radiation. One typical application would be to focus light. Although an electron is considered to be a particle, it also has properties of a wave that can be bent and focused in a manner similar to electromagnetic radiation.

Lens for electrons are not made of glass, but rather use electrical conductors with voltages and currents to create electric and magnetic fields. Some of the ways an electron beam can be controlled include focus, zoom, deflection and rastering.

For example, wire can be coiled into a circular arrangement, and with current flowing through

the wire core, a magnetic field is created that will focus divergent electrons that pass through it into a more collimated beam.

Conductive plates, diametrically opposed across the electron beam path with a potential difference between them creates electric fields that deflects the beam trajectory in a controlled manner.

A typical electron microscope lens system would be composed of a series of many electron optic devices that create both magnetic and electric fields across the electron beam path. As you can imagine, several challenges are present in these systems to optimize the beam alignment and other properties.

For example, the source of the electrons is not always aligned to the “longitudinal optical axis” of the system. Furthermore, the electron travel may not be parallel to the long axis of the system. To solve this problem, most systems incorporate magnetic deflectors to center and align the incoming electron beam. Additional magnetic coils, often referred to as stigmators, are used to help correct for asymmetries in the system. Finally, apertures are placed along the optical axis to block any electrons that are not aligned close enough to the intended beam axis.

### Supporting Systems

#### Physical Structure

Typically, all materials used in constructing the electron and ion guns are compatible with ultra-high vacuum (UHV) technology, which typically utilize pressures even lower than about  $10^{-10}$  torr. It is important that any materials used in vacuum be UHV-compatible and not out-gas (release molecules when put into ultra-high vacuum).

In addition, internal surfaces are typically smooth to not retain but rather facilitate evacuation of atmospheric gases and water vapor, with this material also non-volatile and non-porous to minimize leaching of gases. Materials that are bakeable (i.e., heated to 150 -

Vacuum Range	Pressure (mmHg, torr)	Molecules/cm <sup>3</sup>	Mean free path
Ambient pressure	759.8 (~1.0 atm)	$2.7 \times 10^{19}$	68 nm
Medium vacuum	$8 \times 10^{-1}$ - $8 \times 10^{-4}$	$10^{16}$ - $10^{13}$	100 mm
High vacuum	$8 \times 10^{-4}$ - $8 \times 10^{-8}$	$10^{13}$ - $10^9$	10 cm - 1 km
Ultra-high vacuum	$8 \times 10^{-8}$ - $8 \times 10^{-13}$	$10^9$ - $10^4$	1 - $10^5$ km

Notes: torr x 133 = Pascal, torr /51.715 = PSI.  $10^{-9}$  torr ~  $10^{-7}$  pascal,  $10^{-7}$  torr ~  $10^{-5}$  pascal  
 Data adapted from: <https://en.wikipedia.org/wiki/Vacuum>

Table 1. Examples of vacuum ranges and values. Data adapted from <https://en.wikipedia.org/wiki/Vacuum>

350°C) help to further liberate gases and enhance/enable high vacuum.

To enable these high vacuum conditions, the gaskets used between connected vacuum flanges need to have low or no volatility and tolerate elevated temperatures. Viton (a synthetic fluorocarbon polymer) or copper are typically used. A ConFlat™ flange (CF) system will typically use a copper gasket ring that is positioned between the round flange surfaces of adjacent components, with each of those flange surfaces also including a circular concentric knife-edge. As the bolts between the flanges are tightened, the knife-edge deforms the copper surfaces to create a seal on both sides of the copper ring.

Feedthrough insulators are made of ceramic. In general, common materials (such as most types of insulated wire, rubber tubing, electrical tape, tape labels, duct tape or anything with adhesive) are not UHV vacuum compatible and must be avoided.

A partial list of UHV compatible materials with very low outgassing rates (torr liter/sec cm<sup>2</sup>) for electron gun components include 1) metals; aluminum and alloys (6061, 2024), titanium, copper (OFHC), silver, stainless steel (304, 316), and tungsten, and 2) ceramics; glazed

ceramics such as porcelain, and glasses such as Pyrex.

Other considerations in the design of UHV chambers include highly polished surfaces, minimally contoured interiors, and minimal welds, with geometry optimized for maximum internal access and a usable volume with minimal optimized surface area that allows a high density of experimental apparatus.

Kimball Physics has a product line of vacuum chambers with multiport CF interfaces that are fabricated with precision geometries and generally milled from monolithic pieces of either 316L stainless steel or various grades of titanium. Both the internal and external surfaces are highly polished. This system includes standardized CF mounting interfaces and incorporates mounting features: 1) external channels and bolt holes for supporting the system and various attachments, and 2) internal grooves circumferentially at most inner port diameters for supporting and positioning the experimental components. Kimball Physics also has a complete line of eV parts, essentially a collection of UHV compatible metal and ceramic parts (an “erector set” for scientists) that can provide structure and connectors to create and allow stable and precision support of systems placed within the vacuum chambers.



---

### Vacuum Systems Overview

The quality of the vacuum affects the performance of the electron gun. To create an optimal electron beam, we need to get the air and water gas molecules out of the way.

Typical vacuum for electron beams is in the range of  $10^{-6}$  to  $10^{-7}$  torr (for reference, 1 atmosphere of pressure is  $\sim 760$  torr). For ultra-high vacuum (UHV) applications, at least  $10^{-9}$  torr is needed. Table 1 provides a nice comparison of the molecules in a cubic volume ( $\text{cm}^3$ ) and the mean free path (the average distance traveled by a moving particle between successive impacts or collisions) at atmospheric pressure and different levels of vacuum at room temperature.

The high vacuum (or similarly low pressure) is usually achieved using two types of vacuum pumps: a rough pump and a turbo-molecular pump.

An ion gauge, which is a miniature ion gun and detector system in itself, is used to measure the gas molecules present in the vacuum chamber and monitor the chamber pressure. The ion gauge is typically positioned in the vacuum chamber near the electron gun, but out of the way of the electron beam and target.

It is essential to establish a stable vacuum. An electron gun with a standard refractory metal cathode requires a vacuum of  $10^{-5}$  torr or better, while a gun with a BaO (barium oxide) or a LaB<sub>6</sub> (lanthanum hexaboride) cathode requires  $10^{-7}$  torr. With some materials, such as Y<sub>2</sub>O<sub>3</sub> (Yttria coated Iridium) cathode, a lower vacuum of only  $10^{-4}$  torr is acceptable.

A poor vacuum can damage the cathode or reduce its lifetime. For example, the presence of oxygen or water vapor while the refractory metal cathode is hot (over 700 K) will cause oxidation of the metals, while at higher temperatures (over 1200 K) nitrogen will further degrade the cathode. Thus, the vacuum must be well established before the gun is turned on, and after the gun is turned off, the cathode must be allowed to cool down (usually for at least a half hour) before the vacuum system is vented.

If the electron gun is energized without a proper vacuum, it can be damaged by arc-overs.

Maintaining a suitable vacuum state also keeps the cathode electron source surface from oxidizing and altering its emitting characteristics.

### Power Systems Overview

Power supplies at Kimball Physics typically feature a modular design with miniaturized power supply clusters, optically isolated signals, and a digital interface controller. A LabVIEW™ software virtual control interface is optionally available for many of the systems. Beam stability is an important requirement with stable energy and emission current controls from the power supply required to achieve this.

Figure 4 provides a schematic block diagram of the power supplies (represented as circles) that apply voltages to the parts of the gun to produce and control the electron beam in this one electron gun example.

The power supply provides multiple voltage and/or current sources to the various elements of the electron gun system. For example, a separate power supply module provides for Beam Energy, Source (ECC-Emission Control Current), Grid (Wehnelt), 1st Anode, Focus and X and Y deflection are present. Typical ranges for each component power supply requirements are shown in the table below the schematic (figure 4).


### Detectors

Monitoring the output of the electron gun is typically used during initial gun operation to position the spot, or align the gun, and to determine various power supply settings for optimum performance. The electron beam can be visually monitored by observing the resultant spot on a phosphor screen or can be electrically monitored by measuring the beam current at the target (Faraday cup). The most desirable arrangement is to have both visual and electrical monitoring of the electron beam.

Some of the possible monitoring choices include:

- A Faraday cup and grounded phosphor screen, mounted side by side, on a linear manipulator. Either the phosphor screen or the Faraday cup is positioned in the beam line using the manipulator.

- A grounded phosphor screen with a centrally located hole and a Faraday cup mounted directly behind the hole. Deflection is used to position the spot on the screen when visual monitoring is desired. When electrical monitoring of beam current is desired, the beam is deflected to aim it through the hole into the Faraday cup.
- An electrically isolated phosphor screen connected to a pico-ammeter that enables view and measurement of the beam current directly off the phosphor screen.

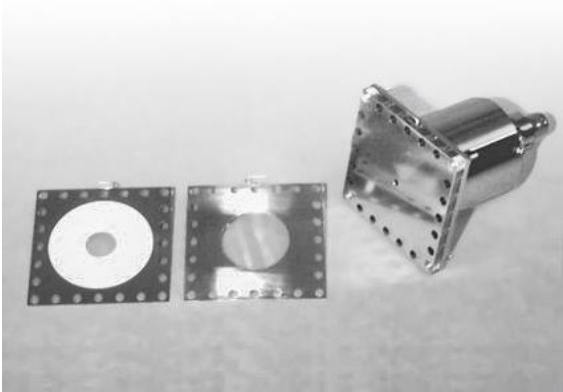

**Figure 4.** Faraday cup with optional phosphor screen.

## Electron and Ion Gun Systems (201)

This section will provide more content about the electron/ion beam and its associated systems to enable the reader to better understand the technical details provided in the technical documents and specifications available for the various electron gun systems.

### Beam Energy

Typical ranges for electron beam energies span several orders of magnitude from 1 eV to 100 keV (100,000 eV). One of the most important characteristics of the beam is the kinetic energy of the electrons or ions. The electron beam energy is determined by the overall potential difference and distance between the cathode and anode that accelerates the electrons as they travel from the cathode source to the final


aperture of the electron gun. This energy, measured in electron volts (eV or keV) and is controlled by the “Energy” power supply channel and can range from 1 eV to 100 keV depending on the gun. One eV is approximately  $1.602 \times 10^{-19}$  joules and is typically the amount of energy 1 electron acquires by accelerating from rest through a potential difference of about 1 volt.

The *First Anode (G-2)*, the first of possibly multiple anodes in the beam path, is referenced several hundred to thousands of volts more positive than the cathode. The electric field created by the potential difference between the cathode and the anode also influences the degree of electron extraction from the cathode. This effect is modified by the electric field present between the Wehnelt or grid and the cathode, which can also modify the number of electrons emitted from the cathode.

The energy spread of the beam is a measure of the variation in energies (typically less than half an eV), as not all particles will exit with the same kinetic energy. The energy spread can vary with different cathode sources. For example, with a 100 keV system, the energy spread for a tantalum (Ta) cathode is ~0.5 eV, Yttrium oxide cathode ( $Y_2O_3$ ) is ~0.4 eV, and Lanthanum Hexaboride ( $LaB_6$ ) cathode is ~0.4 eV.

### Emission and Beam Current

Where the electron goes once emitted largely depends on its energy and the electrical fields in the space around the cathode. If the electron does not have enough energy, or if the space around it has too much negative charge (space-charge limited), the electron will return to the cathode. The electrons that do leave the cathode can be focused and accelerated into the electron beam.

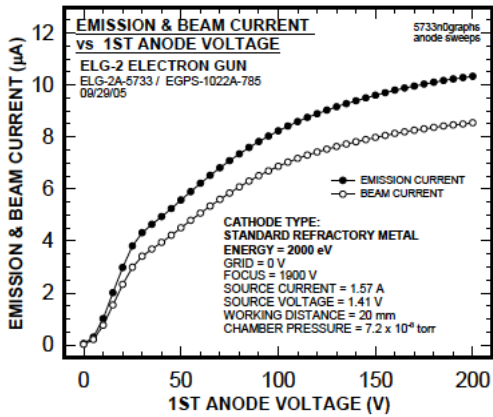



**Figure 6.** The effect of the 1st Anode Voltage on Emission Current from the cathode and the final Beam current.

The emission current is all the electron current that leaves the cathode and eventually goes to ground. This *emission current* depends on both the cathode temperature and the electron beam energy set by the Energy power supply channel (approximately the potential difference between cathode and anode). However, not all the electrons emitted from the cathode will be in the electron beam that finally exits the gun (this can be measured by a Faraday cup at the end of the gun).

Some of the electrons may land on other elements in the gun column and return to ground, but not through the electron beam. Thus, the final electron *beam current* at the target will be less than the initial *emission current*.

As shown in electron gun example Figure 6, increasing first anode voltage increases the emission current, and thus increases the final beam current for a given set of operating parameters. This effect levels off at higher anode values, reflecting the change from a space charge limited situation to a temperature limited one. The first anode potential does not determine the final energy of the electrons, which is reached in this example, only at the grounded second anode after passing through the focus region.

The Wehnelt or grid voltage (G-1) will also affect the beam current. Increasing the Wehnelt potential makes the Wehnelt aperture more negative with respect to the cathode. As the Wehnelt potential increases (becomes more negative, also called increasing the grid bias), the electric field between the cathode and Wehnelt suppresses electron emission from the cathode perimeter, leaving only the center of the

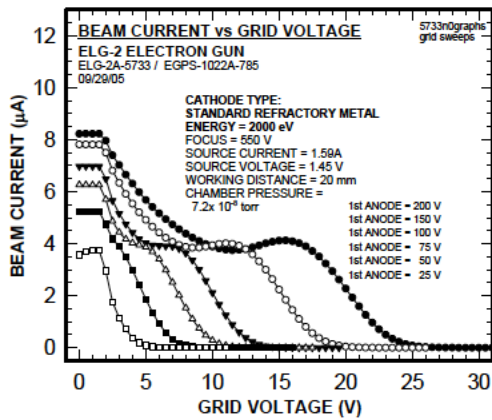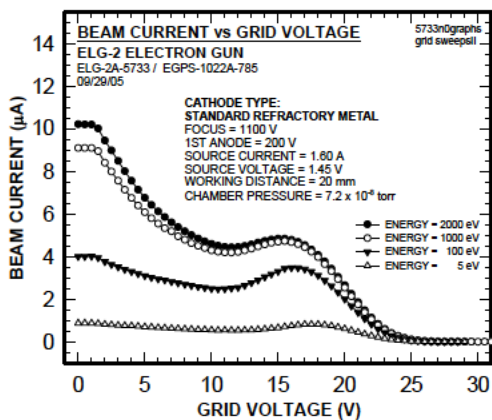

**Figure 7.** The effect of Grid Voltage on final Beam Current, showing cutoff at different Energy levels on the left, and showing cutoff at different 1st Anode voltages on the right.

---

cathode to emit. If enough Wehnelt potential is applied, the beam will be completely suppressed, and the beam is said to be cut off.

This cut-off effect of increasing grid voltage is illustrated in Figure 7 at different Energy levels. Although Energy level affects the amount of beam current, it does not affect the cut-off point on this example system. The figure also shows the effect of different first anode (G-2) levels on beam suppression by the grid (G-1); the grid cutoff value is an increasing function of anode voltage.

Grid cut-off can be used to pulse the electron beam off and on. The various ways that pulsing can be accomplished using the grid are described in the pulsing sections below.

The electric field created by the Wehnelt also controls beam divergence and uniformity by varying electron trajectory.

#### Beam Spot Size

The spot size is the diameter (cm, mm or  $\mu\text{m}$ ) of the electron/ion beam at a given distance from the gun (working distance). This may range from a few microns to several hundred millimeters, depending on the gun.

The spot is measured either visually on a phosphor screen, or by a Faraday cup array; as the edge of the circle may not be clear cut, the full-width- half-max measurement (FWHM) is defined as the width that includes all beam current densities greater than half the maximum density.

In a *flood gun*, which does not have focusing lenses, the size of the spot is mainly dependent on the working distance, or how far away the target is from the end of the gun, although other operating parameters can also effect spot size.

#### Working Distance

In setting up the system, the working distance (the distance from the end of the electron gun to the target) will need to be determined. Many of the gun specifications are affected by the working distance; for example, at a longer working distance, there is more effect of noise, and beam current may be less, or the spot size may be larger. In general, the minimum working distance possible for the experiment should be chosen to reduce problems associated with

transport of electrons, including space charge spreading of the beam and beam distortions from stray magnetic fields and stray electric fields (i.e., the Earth's magnetic field and insulator charge-up).

#### Beam Divergence

Beam divergence is the angular spread of the electron/ion beam from the final aperture of the gun (in degrees) at the working distance and ultimately affects the spot size.

#### Beam Uniformity and Current Density

Beam uniformity describes the distribution of beam current within the spot. The beam current can be further characterized by measuring the beam current density- the beam current per unit area at the target (typically in  $\mu\text{A}/\text{cm}^2$ ).

The variation of beam current density at different positions within the spot is called the *beam current distribution*. In most guns, this is a Gaussian distribution; the spot is brightest in the central area and fades off at the edges. Other electron guns have electron optics elements designed to produce a flatter, more uniform electron distribution across the entire spot.

#### Brightness

Brightness or source brightness is generally the electron density or number of electrons per unit area or (spherical sector) or is commonly defined as the ratio between the beam current and the beam emittance. *Emittance* is a property of a charged particle beam and is a measure of the average spread of particle coordinates in position and momentum phase space, typically with dimensions of length or length times angle. Increasing brightness is often achieved by increasing beam current with reduction in transverse emittance.

High brightness applications (which require small spot and high resolution) include scanning electron microscopes, transmission microscopes, electron probes, scanning Auger systems and electron lithography systems.

#### Electron Optics


Below we continue our discussion about the electron optics elements that modify the transverse motion and axial velocity of the electron beam particles. The electron has an

electrostatic charge and therefore its path can be deflected by an electric field. Furthermore, the electrons in the beam are moving, similar to the current in a wire, whereby there is a magnetic field associated with the beam and therefore it can be deflected by an applied magnetic field. The particles typically have minimal velocities perpendicular to their primary direction of motion. Thus, both electrostatic and magnetostatic fields can both be used to provide forces to change the properties and path of the electron beam. This includes beam focusing, zoom, deflection, rastering, stigmators and condensers. Another interesting property of the electron is that they will release energy as they change direction (i.e., release of photons in a free-electron laser).



#### Beam Focusing

Focus controls the diameter of the electron beam as it approaches the grounded aperture at the end of the gun. In a focusable gun, additional elements are present in the gun to further control the size of the beam.

One focusing approach consists of three axially aligned cylinders which together form an electrostatic converging lens called an Einzel or single lens. The outer two cylinders are typically grounded while the middle cylinder is held at a fixed voltage (Focus Power supply). The electrostatic field is symmetric (see Figure 8) so that if the particles will regain their initial energy on exiting the lens although the direction of the outer particles will be changed such that they converge on the axis.

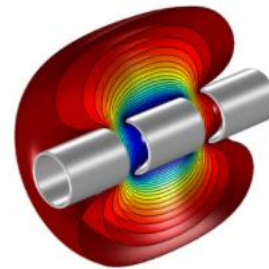

**Figure 8.** Computer model of the electrostatic field of an Einzel Lens for focusing charged particles. [Reference](#)

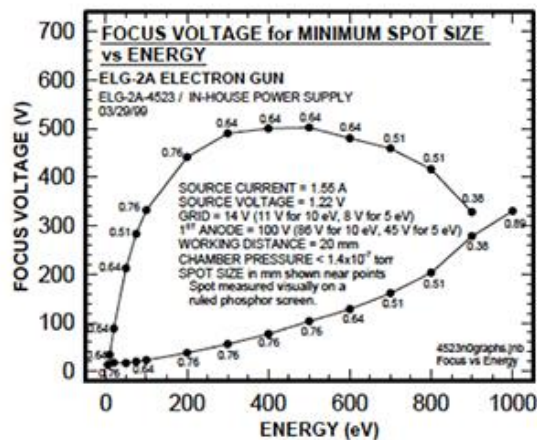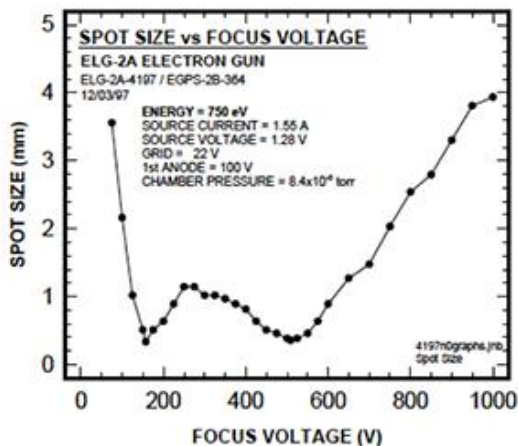

**Figure 9.** The effects of focus on spot size at a particular energy, showing two minimums (left), and The Lens Operating Curve, with the focus values required to produce the minimum spot size (right).

Some guns will have more complex lens systems, such as the zoom lens (see below), for focusing and accelerating or decelerating the electrons. The focus element is controlled by a voltage source referenced to the negative energy supply, typically in the range of 0 to 2 kV. Increasing the focus potential makes the focus element more positive with respect to the cathode. Focus, in conjunction with potentials on the other elements, affects spot size and forms part of the zoom lens. Focus controls the diameter of the electron beam as it approaches the grounded aperture at the end of the gun.


### Zoom Lens

The Zoom Lens consists of the electric fields created by the potentials on first anode elements, focus and the grounded aperture components prior to deflection (please see figures 2 and 4 above). The beam is pre-accelerated in the first anode section and then focused and accelerated or decelerated in the zoom lens section. The beam attains its final kinetic energy at the grounded aperture, the second anode, which completes the zoom lens.

The zoom lens allows flexibility of control because the ratio between the voltages of each element can be varied; in addition, there can be less aberration with more lens elements. Beam energy, beam current, and beam divergence are all directly controllable.

The ratio of Focus voltage to 1st Anode voltage roughly determines the spot size, while the ratio of Energy to 1st Anode voltages determines the amount of acceleration or deceleration that the electron beam undergoes. Since Focus and 1st Anode are referenced to Energy, this lens configuration allows spot size and beam current to remain roughly constant as energy is varied.

The initial elements of this electron gun example all float on the Energy supply; electron energy in this portion of the gun is determined by the potential difference between the cathode and first anode, and as modified by grid and focus. As the electron beam leaves the floating region of the gun and approaches the grounded aperture, it is either accelerated or decelerated, depending on the relative voltage differences between the Energy supply and the 1st Anode supply.

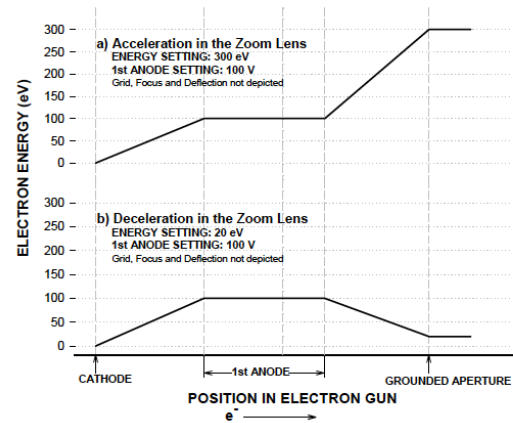


**Figure 10.** Example of Zoom Lens Acceleration / Deceleration Schematic

A theoretical schematic energy profile over the length of the gun might be illustrated by Figure 10. For example, if Energy is set to 300 V and 1st Anode is set to 100 V, then the electron beam encounters potentials from -300 V to -200 V to 0 V (relative to ground potential) and attains energies from 0 eV to 100 eV to 300 eV as it passes through the gun. As illustrated in Figure 10(a), the beam is accelerated from 100 eV to 300 eV in the final zoom lens portion of the gun.

Alternatively in Figure 10(b), if Energy is set to 20 V and 1st Anode is set to 100 V, then the electron beam encounters potentials from -20 V to +80 V to 0 V (relative to ground potential) and attains energies from 0 eV to 100 eV to 20 eV as it passes through the gun. The energy profile for this case illustrates that the beam is decelerated from 100 eV to 20 eV in the final zoom lens portion of the gun. If the deceleration ratio (1st Anode to Energy) exceeds 10 to 1, then spot size and beam current may deteriorate.

### Deflection / Rastering

The deflection elements consist of two pairs (X and Y) of deflection plates located downstream of the focus element. Potentials applied to these plates produce an electrostatic deflecting force in a plane perpendicular to the direction of electron travel, allowing the beam to be aimed or rastered (for units with the Raster Generator option). Deflection allows the beam to be guided by the operator to the target.

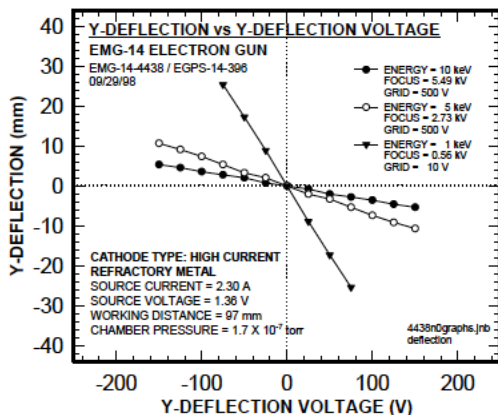

**Figure 11.** The effect of deflection voltage on the position of the spot a three different energy in a typical electron gun example.

The position of the spot varies directly with the deflection voltages applied (as noted in the Figure 11). However, the amount of movement also depends on beam parameters such as the acceleration energy. At low energies, a given deflecting voltage produces a much larger shift in position. The maximum possible deflection may be less at higher energies than at low energy.

Non-uniformities in the deflecting electric field present at the edges of the deflection plates can cause the spot to become distorted if it is deflected too close to the deflection plates. Distortion varies with both the diameter of the electron beam as it passes through the deflection plate region and the potentials applied to the deflection plates. The greater the diameter of the electron beam as it passes through the deflection plates, the greater the distortion.

Electron beam diameter in the deflection plate region can be minimized by choosing a focus voltage which places the beam crossover between the deflection plates. This however will not minimize spot size at the target.

If a minimum spot size is desired, then a focus voltage must be chosen which places the beam crossover at the target. Under these conditions, the beam diameter as it passes through the deflection plate region is not at a minimum and

there will exist more deflection distortion. Better spot shapes will be achieved if the amount of deflection is kept to a minimum.

### Magnetostatic Lens

A magnetic lens typically consists of several electromagnets (essentially coils of wire) placed at the vertices of a polygon in a plane (perpendicular to the axis of the particle beam) to create a customized magnetic field to manipulate the particle beam. For example, four, six and eight electromagnets would create quadrupole, hexapole and octupole magnetic lens systems.

Magnetic beam rocking, which is used in a few high-energy guns, is very similar to rastering. Instead of electrostatic deflection plates, currents in electromagnetic coils around the gun move the electron beam in an unsynchronized pattern over the target area. The effect is the same as rastering, to provide a more uniform coverage of a larger area.

The passing particle is subjected to two vector forces, with one parallel to the core and the other parallel to the lens radius. The magnetic field parallel to the lens radius causes the particle to spiral which then allows the field parallel to the core to influence and focus the electron. The magnetic fields are typically not homogeneous, with particles close to the center less effected than the particles farther away from the beam axis.

### Cross Over Points

Along the beam path there are typically several cross over points. A cross over point is simply a location in the electron beam path where the converged beam reaches a minimum size or cross-section. The brightness or beam current density will typically be greatest at this point. The first cross over point is typically just past the Wehnelt cylinder.

### Beam Operation and Performance Factors

The section below speaks about the temporal control of the particle beam in including methods

---

to turn the beam on and off using the Wehnelt or grid potential along with the various methods available to pulse the beam.

#### Beam On/Off

In most guns, the electron or ion beam may be turned off and on while the gun is running. The way this is accomplished depends on the specific gun design; often several methods are available for an electron gun and some of them are discussed below.

#### Grid Cut Off

The grid, sometimes called G-1 or the Wehnelt, is the first electrostatic element in the gun that the electrons/ions encounter after being emitted from the cathode/ion source. The grid serves as the first control point over the beam and usually can be used to shut off the beam.

In an electron gun, if the grid voltage is sufficiently negative with respect to the cathode, it will suppress the emission of the electrons, first from the edge of the cathode and at higher (more negative) voltages from the entire cathode surface. Please see Figure 7 above for an example of Grid Voltage vs Emission Current.

Similarly, in some ion guns, a positive grid voltage can suppress the positive ions from the ion source. The minimum voltage required to completely shut off the flow of electrons/ions to the target is called the *grid cut off*. The grid voltage is typically manually controlled by a potentiometer knob on the power supply; thus, in most guns, the beam can be turned off while the gun is running by setting the grid to the “cut off” voltage. Other automated methods to control the grid voltage are covered below.

#### Beam Pulsing

Pulsing, also called fast beam pulsing, is stopping and starting the flow of electrons or ions in a fast cycle. This pulsing is usually accomplished by rapidly switching the grid voltage to its cut off potential to stop the beam. The grid voltage can be controlled by several different methods (listed in order of speed):

- Manual control with the dial potentiometer
- Remote control with computer input into power supply terminals

- Dual grid power supplies controlled with a TTL signal input
- Capacitive pulsing with a pulse junction box or cable and an external pulse generator.

The simplest method of turning the beam off and on is just to cycle the grid voltage by hand with the control knob on the front of the power supply. Clearly, this would be slow and not reproducible.

A more systematic method of controlling the grid is by an input signal into the remote terminals on the rear panel of the power supply. Remote control is a standard feature on all our Kimball Physics power supplies, so this method does not require any system options. However, it may not provide sufficiently fast pulsing.

With the dual grid pulsing option, there are two grid power supplies built into the main power supply. A pulsing TTL (transistor-transistor-logic) signal switches rapidly between the two supplies, pulsing the beam on and off. For most guns, the dual supplies are (1) the normal, variable control grid supply which is adjusted to allow the electron flow and (2) a fixed grid supply which is fixed at the cut-off grid voltage at the factory.

For guns that usually have a positive grid, the dual supplies are (1) a variable positive grid supply which allows the electron flow, and (2) a variable negative grid supply which is adjusted to cut-off.

For the capacitive or fast pulsing option, many guns can be equipped with a capacitor-containing device (either a separate pulse junction box or a cable with a box) that receives signal from an external pulse generator. The grid power supply and pulse generator outputs are superimposed to produce the voltage at the grid aperture. The grid power supply is first set at the cut off voltage. As the grid voltage is negative with respect to the cathode, positive voltage pulses are required to pulse the gun on. The general pattern of the beam pulsing is a square wave with a variable width (time off and time on) and a variable repetition rate. Capacitive pulsing can provide the fastest rise/ fall time and shortest pulse length of the various methods. However, the capacitor does not permit long pulses or DC operation.



---

## Electron Sources

Electron guns/sources are often classified by the type of electric field generation (DC or RF), the emission mechanism (thermionic, photoelectron, cold or field emission and plasma source), and by the number of source electrodes. Our discussions will focus on DC field generation, thermionic emission, and single cathode systems.

As previously introduced, the simplest source of electrons in an electron gun cathode could be a metallic wire or filament such as tungsten bent into a hairpin shape, where we apply a small voltage to pass a current through the wire's resistance to heat the wire. However, most electron emitter configurations used by Kimball physics indirectly heat the cathode material by thermal conduction from the surrounding support material. The support structure is heated by placing a potential difference (~ 1 - 4 volts) across its resistance, with a "source" current (~1.5 - 3A) heating the conductor.

Kimball Physics emitters provide a reliable source of electrons or ions for use in a variety of instruments including scanning and

transmission electron microscopy, electron lithography, x-ray generation, free electron lasers, electron accelerators, and other custom applications. Other applications include our own Kimball Physics electron guns and ion guns.


Cathode types for electron emission available at Kimball Physics include:

- Lanthanum Hexaboride (LaB<sub>6</sub>) high brightness single-crystal
- Tantalum (Ta) disc
- Barium Oxide (BaO) coated disc
- Ytria (Y<sub>2</sub>O<sub>3</sub>) coated iridium disc

Cathodes are available on several types of bases to interface with multiple systems:

- Standard ceramic AEI base
- Compact Kimball Physics CB-104 base
- Compact Kimball Physics CB-105 base
- Various microscope bases
- Cathodes may also be mounted on customer-supplied based

Please refer to our cathode documents for more thorough discussions of Kimball Physics various cathode parameters and products.

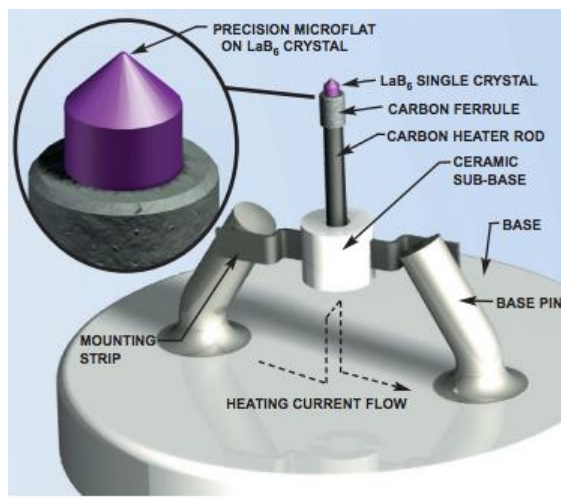

**Figure 12.** Lanthanum Hexaboride (LaB<sub>6</sub>) cathode model with heating current path through slotted precision machined single-piece carbon rod and mounting strips with sub-base for rigidity and easier mounting. The diameter of the LaB<sub>6</sub> prior to the conic section is about 440 microns. Cone angles are typically 60 or 90 degrees with a microflat at the tip ranging from 6-300 microns.

---

## Other Types of Electron Emission

### Photoelectron Emission

Also, known as the photoelectric effect, it characterizes the emission of electrons from a material surface when light falls on that object. It involves the transfer of energy from the light to the electron to induce changes in the kinetic energy of the electron.

The energy in the photon to dislodge the electrons must reach or exceed a threshold energy. Below this threshold, no electron will be emitted, even with an increase in light intensity (number of photons per unit time) or exposure time. For example, emission of conduction electrons from typical metals requires a few electron volts (eV), corresponding to short wavelength (high frequency) or ultraviolet light where the photon energy ( $E = hv$ ) is related to the product of Planck's constant  $h$  and the light frequency ( $\nu$ ).

This photoelectric effect can occur in any material. However, as electrons are emitted and a charge imbalance is created, the conductivity of the material determines how much current (electrons) can flow into the material to neutralize the charge imbalance. The energy barrier to photoemission is usually also increased by thin oxide layers on the material surface. Metals with clean surfaces and minimal exposure to oxygen in a vacuum environment are more favorable emitters.

### Field Emission

Field emission, also known as field electron emission, is the emission of free electrons induced by a strong electrostatic field. The most common configuration would be from a sharply pointed solid surface into a surrounding high vacuum. Field emission in pure metals requires

high electric fields ( $\sim 1$  gigavolt/ meter), high vacuum ( $\sim 10^{-10}$  torr) and are also strongly dependent on the work function of the material.

Applications that would benefit from high brightness electron sources include high resolution microscope and charge neutralization applications (spacecraft). Cold field electron emission usually involves minimal heating of the cathode, with an initial state is internal thermodynamic equilibrium. The geometry of the field emitter tip is typically a sharply pointed wire that enables smaller spot sizes and higher current densities. Methods are also available to thermally assist with field emission and lower the magnitude of the electric field required.

## Ion Sources and Systems

There are three basic processes by which ions are generated in the several Kimball Physics ion guns: electron impact gas ionization, microwave gas ionization, and alkali metal solid surface ionization.

### Gas Ionization

This type of ion source has a filament which emits electrons when heated by the Source power supply. An inert or a reactive gas, such as argon or oxygen, is introduced from an external tank via a gas feedthrough into the region inside the ion gun near the filament. The electrons emitted from the filament are accelerated into the gas region and collide with the neutral gas molecules. If the impact is of sufficient energy, an electron will be removed from the neutral gas molecule, resulting in a positive ion. Usually, the ion will be singly charged. Depending on the energies involved, the primary and the secondary (released) electrons may go on to collide with other gas molecules and cause further ionization. The electrons finally land on a positive structure called the ion cage.

---

The number of ions produced by electron impact ionization depends mainly on the number of electrons emitted, their energy, the type of gas and the number of gas molecules present to be ionized. The ionization potential of each gas species determines the minimum energy the bombarding electrons must have to remove an outer electron. Within a normal range, the pressure of the gas introduced into the gun is linearly related to the ion beam current produced.

#### Plasma

The microwave ion source is similar to the electron impact ion source in that it generates ions from a gas introduced into the gun. However, instead of electron bombardment, it uses microwave power, which is transmitted from an external generator to a helical antenna in the ionization chamber, to heat the gas and create a plasma, a highly-energized state in which the gas molecules are disassociated into ions and free electrons. The plasma can be seen as a glowing gas in the chamber, similar to a fluorescent light tube. After the ions are extracted by several apertures, they pass through an ExB filter that consists of a region with an electrical field (E) perpendicular to a magnetic field (B). This ExB filter separates the ions based on their mass; by varying the electric field, ions of different masses can be selected as in a mass spectrometer. The electrical field is adjusted so that only ions of the type desired (based on their mass) can pass through, while other ions and molecules are deflected. Further lenses control the energy, focus and deflection of the selected ions.

#### Solid Ionization

In other ion guns, ions of alkali metals such as  $\text{Cs}^+$ ,  $\text{Li}^+$ , or  $\text{K}^+$ , are generated directly by solid surface ionization. Instead of having a filament, this type of ion source has a cartridge containing solid compounds of the alkali metal. When heated by the source voltage, these compounds undergo a solid-solid chemical reaction that releases the alkali metal ions at the surface.

These positive ions are then accelerated and focused into a beam as in other guns.

#### Ion Beam

The ions generated by the ion source are then formed into the ion beam. A gun element called the extract aperture, which has a negative voltage relative to the ion energy, accelerates and extracts the positive ions from the region of electron bombardment; this is similar to the function of the anode in the electron gun triode. The grid element, with its potential controlled by the Grid power supply, also helps extract and focus the ions. Depending on the gun, additional elements may be present to accelerate, decelerate, focus and deflect the beam.

---

## Frequently Asked Questions- Gun Operations

### HOW IS THE BEAM MOVED ON THE TARGET?

#### Mechanical Alignment

Mechanical alignment is available with many gun mounting systems to aid in aiming the electron or ion beam at the target. A rotating flange or a device consisting of two rotating angled disks (a Port Aligner) can be used to mechanically change the position of the gun. Port Aligners are available from several manufacturers.

#### Deflection

Many Kimball Physics electron and ion guns have an electrostatic deflection system as an option. In the gun column following the focusing lenses, two pairs of plates (X+, X- and Y+, Y-) are arranged around the path of the beam. Voltages are applied to these plates by manually controlled X and Y power supplies, and the resulting electrical field bends the path of the beam. This deflection is used to move the beam around in the plane of the target to position the beam. In a few guns, magnetic coils are optionally available to similarly deflect the beam for alignment.

#### Rastering

Rastering, which is related to deflection, is a continuous movement of the electron or ion beam over the target plane to uniformly cover an area over time. This rastering is similar to what occurs with a TV or CRT screen. Two synchronized, cyclically varying voltages are applied to the X and Y deflection plates in the gun. This causes the beam to move in a synchronized pattern: sweeping back and forth (X direction) while simultaneously moving down more slowly (Y direction), then returning to the start position without sweeping back and forth (retrace). Visually, it appears that the beam covers a large square area, instead of a single small spot. Rastering is occasionally done in an unsynchronized mode.

#### Beam Rocking

Magnetic beam rocking, which is used in a few high-energy guns, is very similar to rastering. Instead of electrostatic deflection plates, currents in electromagnetic coils around the gun move the electron beam in an unsynchronized pattern over the target area. The effect is the same as rastering, to provide a more uniform coverage of a larger area.

### HOW IS THE ELECTRON OR ION BEAM OBSERVED?

Because an electron or ion beam is invisible, it must be observed by its effects, by employing some type of detector. Two common types of detectors are phosphor screens, which emit light when struck by the electrons/ions, and Faraday cups, which collect the electrons/ions and allow current measurement. A variety of phosphor screens and Faraday cups available from Kimball Physics.

#### Phosphor Screen

A grounded phosphor screen in the target position allows visual, real-time observation of the spot produced by the beam. The phosphor screen emits light (photons) when bombarded by high-energy particles, either electrons or ions; the light color, typically blue or green, depends on the phosphor material. On the screen, the electron/ion beam usually appears as a small, solid circle (the spot). If the beam has not been focused, the spot may be much larger, or have fuzzy edges, or be made up of many dots (beamlets), or even (with an LaB<sub>6</sub> cathode) have an unusual cross shape. As the position, size and shape of the spot can be easily observed, the screen is useful for alignment and setting the proper operating parameters for the gun.

The use of ion beams with phosphor screens has been challenging in that they often become contaminated quickly with a shortened lifespan. The use of a Microchannel Plate (array of electron multiplier tubes) in front of the phosphor screen seems to help mitigate this problem.

---

### Faraday Cup

A Faraday cup can be used to detect and measure the actual beam current emitted from the electron/ion gun. The Faraday cup consists of a shielded cup with an aperture, which collects the electrons or ions, and an output wire, which is connected to an ammeter. The Faraday cup can be completely separate from the gun, or it can be part of an assembly mounted to the end of the electron/ion gun and manipulated remotely. Some Faraday cup assemblies include a phosphor screen as well. An array of small Faraday cups or a Faraday cup on mounted on a linear manipulator can be used to measure the distribution of the beam current across the spot; this shows the beam uniformity.

### HOW ARE THE POWER SUPPLIES CONTROLLED?

The power supply unit, which is part of the electron or ion gun system, is a complete unit containing electronic circuitry, meters and controls, that takes power input from the outside line and converts it to voltages and currents required to run the electron or ion gun. Individual power supplies within the power supply unit provide the voltage to specific gun elements. The output of a power supply is monitored by either an analog meter (one with a dial and pointer) or with a digital voltage meter (DVM).

### Manual and Remote Control

These individual supplies can be controlled either manually by the controls on the front of the power supply unit, or remotely by a signal into the remote controls on the rear of the power supply. The manual dial potentiometer controls of the FlexPanel directly controls the output of the individual power supply for the various electron gun elements. By use of the remote-control input terminals, the supplies may be either voltage programmed or resistance programmed. The voltage control signal can be obtained from any source, such as a computer with a digital to analog converter, or an analog supply. Resistance programming involves connecting a resistor to the remote terminals.

### Feedback Control (ECC)

In addition to the usual manual or remote control, the source power supply can be controlled in a

feedback-regulated mode, labeled ECC for Emission Current Control. In electron guns, this feedback circuitry regulates the source power supply, which heats the filament, so that a constant emission current is maintained from the cathode. Similarly, in alkali metal ion guns, the ECC regulates the heating of the ion cartridge and so maintains a constant ion emission current. In gas ion guns, the ECC feedback circuitry controls the electrons used for ionization, but this does not completely regulate the emission of ions from the ion source, as the ionization is also dependent on the gas pressure and other factors.

### HOW IS THE GUN CONNECTED TO THE VACUUM SYSTEM?

#### Mounted Guns

Most Kimball Physics electron and ion guns are designed in a mounted configuration, the gun being housed in a tube with a flange for connecting the vacuum system. The mounting system uses a standard type of vacuum flange (a CF or ConFlat flange) that has a knife edge on the sealing surface and comes in standard sizes according to its approximate diameter; for example, a 1.33" CF flange has about 1.33 inch or 34 mm diameter, a 2.75" CF flange (70 mm), and a 6.0" CF (152 mm).

The mounted gun also includes cable connectors for the power supply and vacuum feedthroughs. These are insulated connections for the electrical wires and other tubes needed to connect the gun to the power supply cable and to other external parts, such as a gas tank.

A Kimball Physics Flange Multiplexer, which is used as the mounting system on many electron/ion guns, has a one (2.75" CF) flange surface to connect to a vacuum port and five smaller (1.33"CF) flanges for various feedthroughs and cable connectors.

#### Unmounted Guns

Some of the simpler guns are also available in an unmounted configuration, which allows the customer to design the housing, mounting system and vacuum connections.

---

#### HOW ARE THE GUNS TESTED?

Every new electron or ion gun and power supply, built at Kimball Physics, is individually tested by being run in our vacuum laboratory, with the data on various gun parameters collected. This ensures that each gun functions properly, is well aligned, and meets its specifications. Graphs are provided to the customer showing the results of this testing to aid in determining the operating conditions for the particular gun. Guns returned for rebuild or repair undergo similar, but less extensive, testing. (The only exceptions to this testing procedure are a few simple, firing unit models sold in large batches.)

#### Data

The data collected depends on the gun model and the needs of the customer. The following graphs are generally included: 1) the V-I Characteristic (Source Current vs Source Voltage), which shows the performance of the particular cathode/ion source; 2) Emission and Beam Current vs Source Current, which shows the overall gun performance, how much beam can be produced, and is used to set the source when running the gun; 3) Emission Current vs Grid Voltage and 4) Beam Current vs Grid Voltage at various Energies, which together

show how the grid can be used to improve or to cut off the beam.

Depending on the electron gun, graphs showing focusing parameters, spot size, beam current density, or other properties may also be provided. Additional custom-designed testing procedures can be performed as required by the customer.

---

Intentionally left blank.

---

## References

### Lanthanum Hexaboride (LaB<sub>6</sub>)

#### Technical Bulletins:

LaB6-01 Technical Bulletin - Cathodes / Emitters- General Guidelines for Operating ES-423E LaB6 Cathodes.

LaB6-02 Technical Bulletin - Cathodes / Emitters: The Relationship Between LaB6 Cathode Life and Gun Vacuum

LaB6-03 Technical Bulletin - Cathodes / Emitters: Emission Drift- Lab6 and Gun Stability

LaB6-04 Technical Bulletin - Cathodes / Emitters: Oxygen Activation of LaB<sub>6</sub> Cathodes- The Double Saturation Effect

LaB6-05 Technical Bulletin - Cathodes / Emitters- Kimball Physics ES-423E LaB6 Cathode, Style 60-06 (60-degree cone angle, 6 micron Flat)

LaB6-06 Technical Bulletin - Cathodes / Emitters- Kimball Physics ES-423E LaB6 Cathode Operating Instructions for Leica / Cambridge Stereoscan Series SEMs

LaB6-07 Technical Bulletin - Cathodes / Emitters- Recovery of Emission from ES-423E LaB6 Cathodes Following a Vacuum Dump.

#### LaB<sub>6</sub> ES423-E

LaB6 Cathode (ES-423E) - Cathodes / Emitters- Extended Life LaB6 Cathode- Description and Specifications

LaB6 Cathode (ES-423E) - Cathodes / Emitters- Extended Life LaB6 Cathode- User Information.

#### LaB<sub>6</sub> ES423-GR

LaB6 Cathode (ES-423GR ) - Cathodes / Emitters- Extended Life LaB6 Cathode- Description and Specifications

ES-423GR Guard Ring LaB6 Cathode User Information

#### LaB<sub>6</sub> ES440

LaB6 Cathode (ES-440) - Cathodes / Emitters- Extended Life LaB6 Cathode- Description and Specifications

ES-423GR Guard Ring LaB6 Cathode User Information

#### LaB6 Other Information

The Relationship Between LaB6 Cathode Life and Gun Vacuum.

LaB6 Surface Recession vs Temperature - Cathodes / Emitters (Chart/Graph)

LaB6 Surface Recession vs Theoretical Brightness - Cathodes / Emitters (Chart/Graph)

#### Barium Oxide Cathodes

Barium Oxide Cathodes / Emitters- Description and Specifications

Barium Oxide Cathode (ES-015) - Cathodes / Emitters- Care and Operating Instructions

Barium Oxide Cathode (ES-064) - Cathodes / Emitters: Care and Operating Instructions

Barium Oxide Cathode (ES-066) - Cathodes / Emitters: Care and Operating Instructions

#### Tantalum Cathodes

Tantalum (Ta) Cathodes / Emitters- Description and Specifications

Ta Disc Cathode (ES-042) - Cathodes / Emitters: Care and Operation Instructions

Ta Disc Cathode (ES-044) - Cathodes / Emitters: Care and Operating Instructions



---

Ta Disc Cathode (ES-046) - Cathodes /  
Emitters- Care and Operating Instruction

### **Yttria Coated Cathodes**

Yttria Coated Iridium Cathodes / Emitters-  
Description and Specifications

Yttria Coated Cathode (ES-529) Y2O3 Disc  
Cathode Care and Operating Instructions

Yttria Coated Cathode (ES-535) - Cathodes /  
Emitters- Care and Operating Instructions

Yttria Coated Cathode (ES-535W) - Cathodes /  
Emitters- Care and Operating Instructions

Yttria Coated Cathode (ES-526) - Cathodes /  
Emitters- Care and Operating Instructions

### **Tungsten Cathodes**

Tungsten Filaments- Hairpin Thermionic  
Emitters

#### Other References

Hongliang Liu, Xin Zhang, Shuyu Ning, Yixin  
Xiao, Jiuxing Zhang. The electronic structure  
and work functions of single crystal LaB6 typical  
crystal surfaces. Vacuum, Vol 143, September  
2017, Pages 245-250

---

## Glossary

### **AEI Base**

An industry standard base for cathode emitters. Typically, the ceramic disk dimensions are approximately 12.0 mm in diameter with a thickness of 2 mm. The spacing between the center of the pins on the socket side is about 6.45 mm, with 1.0 mm pin diameters and 4.7 mm pin lengths.

### **Alignment**

- 1) Positioning of the physical elements of the electron/ion gun so that the charged particles travel in the optimal path through the gun.
- 2) Mechanical beam alignment: A mechanical adjustment of the gun that changes the orientation of elements of the gun with respect to other elements or with respect to the target; use of a device to line up gun with the target, for example, using a Port Aligner which consists of two rotating angled disks.
- 3) Magnetic compensation alignment: Displacement of the electron/ion beam in a plane (X-Y) perpendicular to the line of travel (Z) of the beam, by means of the magnetic fields produced by electric coils (X+,X- and Y+,Y-) arranged around the beam; used to position the beam on the target and improve beam current and beam uniformity.

### **Analog Meter:**

A meter with a dial and pointer which displays current or voltage.

### **Amperes, Amps; A:**

A measure of current (electron flow).

1 Ampere = 1 Coulomb/sec =  $6.242 \times 10^{18}$  electrons/sec

### **Anode: (also G-2)**

Often referred to the more positively charged electrode

The second electrostatic lens element in an electron gun, which usually provides the more positive potential difference with respect to the

cathode electron source that accelerates the electrons

The third element of the triode (cathode, grid and anode); usually a tube with an aperture (hole); often grounded.

### **Aperture:**

A small hole that the electron/ion beam passes through, which can limit the divergent trajectory of the beam path and also control the diameter of the beam.

### **Auto-bias:**

Most electron guns in SEM and TEM use auto-bias operation, whereby the only electrical variable other than the cathode heating is the value of the auto-bias resistor.

### **Barium Oxide (BaO):**

A coating applied to a cathode electron source that reduces the work function enabling electron emission at lower temperatures, thus also reducing light output.

### **Barium Oxide (BaO) Cathode:**

A Kimball Physics thermionic low light and low temperature electron emitter, with these characteristics related to its lower work function.

### **Bake out:**

Heating the electron/ion gun while in vacuum to remove and pump away contaminants.

### **Beam Blanking:**

Deflecting the electron/ion beam to one side of the gun tube to interrupt the flow of particles to the target without actually turning off the beam; can be used to pulse the final beam current repeatedly on and off, controlled by a TTL signal input.

### **Beam Current:**

The electron/ion current exiting the gun, usually, measured by a Faraday cup at the end of the gun (in mA,  $\mu$ A, or nA); not identical to emission current.

**Beam Current Density:**

The beam current per unit area reaching the target (typically in  $\mu\text{A}/\text{cm}^2$ ).

**Beam Divergence:**

The angular spread of the electron/ion beam as it exits the gun (in degrees  $^\circ$ ).

**Beam Uniformity:**

A qualitative statement about the beam current distribution, usually Gaussian-like.

**Beam Spot Size:**

The spot size is the diameter of the electron/ion beam at a given distance from the gun (working distance). Typically range from 10  $\mu\text{m}$  to several hundred  $\mu\text{m}$ , depending on the gun. The spot is measured either visually on a phosphor screen, or by a Faraday cup array; as the edge of the circle may not be clear cut, the full-width- half-max measurement (FWHM) is defined as the width that includes all beam current densities greater than half the maximum density. In a flood gun, which does not have focusing lenses, the size of the spot is mainly dependent on the working distance, on how far away the target is from the end of the gun.

**Bias Resistor, controllable:**

A resistor used to control emitter heating current more accurately.

**Brightness:  $\beta$ :**

Brightness or source brightness; electron density or number of electrons per unit area or (spherical sector). High brightness applications (which require small spot, high resolution) include scanning electron microscopes, transmission microscopes, electron probes, scanning Auger systems and electron lithography systems.

$$\beta = \sigma / \pi k T$$

$$\sigma = \text{work function} = eV$$

$$k = \text{Boltzmann constant} = 8.617 \times 10^{-5} \text{ eV/K}$$

$$T = \text{absolute temperature} = \text{K}$$

$$\text{Units} = \text{Current/unit area/unit solid angle}$$

**Brightness Limited:**

Brightness-limited electron optical systems: SEM's, TEM's, probes, electron lithography systems, etc.

**Cartridge:**

(1) Firing unit cartridge: For electron Guns: An assembly containing cathode and grid which is part of a larger firing unit of some guns and can be replaced separately.

(2) Ion source cartridge or alkali metal cartridge: For Ion Guns: A cup containing a solid that is heated to produce alkali metal ions directly by a solid-solid chemical reaction, part of the ion source firing unit.

**Cathode:**

The structure in the electron gun that emits electrons, also known as emitter, electron source; the first element of the triode (cathode, grid and anode); simple configuration may consist of a sharply bent wire (hairpin) or a heated wire and an emitting disc; several other types of materials and configurations to optimize performance and live-time parameters.

**Cathode Loading:**

$\text{A}/\text{cm}^2$ ; Typical ranges of 3.8 to 80

**Cathode Poisoning:**

Usually refers to an unintended chemical process involving the surface of the cathode that unfavorable alters its emission characteristics.

**CB-104 base:**

A compact Kimball Physics ceramic cathode base

**CB-105 base:**

A compact Kimball Physics ceramic cathode base.

**CF, ConFlat:**

A type of vacuum flange with a knife edge on the sealing surface; sized by the approximate outside diameter: 1.33 CF ~1.33 inch ~34 mm,

---

2.75" CF ~2.75 inch ~70 mm, 4.5" CF ~4.5 inch ~114 mm.

(The ConFlat flange was invented by Bill Wheeler at Varian; ConFlat is a registered trademark of Varian Associates.)

**Current Density:  $j$ ; (emitted current density)**

Derived for thermionic emission with Richardson-Dushman equation (related to work function and temperature). Amperes per meter<sup>2</sup>, A/m<sup>2</sup>

**Cut off (also Grid cut off):**

(1) Shutting off the flow of electrons/ions to the target; for electrons, by making the grid voltage sufficiently negative that the electrons from the cathode are repelled and do not go through the grid, or similarly for positive ions, by making the grid sufficiently positive that the ions in the source region are repelled.

(2) The minimum grid voltage that reduces the beam current to zero.

**DAQ:**

Data acquisition

(1) Communication between a computer program, such as a LabVIEW™ executable program, and a physical system, such as an electron or ion gun power supply

(2) Also refers to computer boards and cables used to transfer data and control signals to and from the equipment

**Deflection, X and Y Deflection:**

(1) Electrostatic or electromagnetic displacement of the electron/ion beam in a plane (X-Y) perpendicular to the line of travel (Z) of the beam; a) Electrostatic by means of voltages applied to two pairs of plates (X+,X- and Y+,Y-), or b) Magnetostatic by currents applied to deflection coils, arranged around the beam; used to position the beam on the target, generally without affecting the focus and energy

of the beam. Also sometimes called centering or X-Y centering.

(2) The two power supplies that apply voltage to the deflection plates. Also the manual control on the power supply panel that control the deflection power supplies.

**Differential Pumping:**

Maintaining a different vacuum pressure in the electron or ion gun and the main vacuum chamber by connecting separate vacuum pumps to the two areas, usually using a Compact Tee differential pumping connector on the gun.

**ECC:**

(1) Emission Current Control for Electron Guns: Feedback circuitry that regulates the source power supply that heats the filament; used to maintain a constant emission current from the cathode.

(2) Emission Current Control For Alkali metal Ion Guns: Feedback circuitry that regulates the source power supply that heats the alkali metal cartridge; used to maintain a constant ion emission current from the ion source.

(3) Electron Current Control for Gas Ion Guns: Feedback circuitry that regulates the source power supply that heats the filament in the ion source, so that a constant flow of electrons is given off; only indirectly controls the production of ions by the ion source.

(4) Also the power supply panel controls and switch that control the ECC circuitry.

**Electric Field, E:**

Typically, a field around an electric charge that exerts a force on other charges, with attraction between dis-similar charges, and repulsion between like charges. Electric fields are created by electric charges (Gauss's law) or by time varying magnetic fields (Faraday's law of induction). For example, an electric field is

present between the conduction plates of capacitor while maintain a voltage (potential difference) between the plates. With an electron gun, an electric field is created between the source cathode and the anode to accelerate the thermionic emitted electrons into an electron beam. An electric field is also typically created between the Wehnelt and the cathode source to modify or temporarily suppress emission.

-on an atomic scale, an electric field is responsible for the attractive force between the electrons and the atomic nucleus.

-SI units, Newtons per Coulomb, N/C

-Volts/meter, V/m

#### Electrode:

An electrode is typically an electrical conductor used to contact a non-metallic part of a circuit. In an electron gun system, the negative electrode is the cathode, the positive electrode is the anode, and the non-metallic medium is the partial vacuum space. In an EKG system to measure electrical cardiac activity, electrodes are placed opposing sides of the chest to analyze the cardiac rhythm.

#### electron Volt (eV):

• a unit of energy equivalent to  $1.6 \times 10^{-19}$  joules. It is amount of energy gained or lost by the charge of a single electron moving across an electric potential difference of 1 volt.

• 1 eV is equivalent to the energy obtained by a particle with unit electronic charge of ( $1.6021 \times 10^{-19}$  coulombs) accelerated by a 1-volt potential difference and is equal to  $1.6021 \times 10^{-19}$  joules.

- $eV = \text{kg m}^2/\text{sec}^2$

- $eV = 1.602 \times 10^{-19}$  joules

- $1 \text{ eV} = E_{(eV)} = V_{(V)} \times Q_{(e)} = (1 \text{ V}) \times (1.6 \times 10^{-19} \text{ C}) \text{ per electron} = 1.6 \times 10^{-19} \text{ J}$

- $1 \text{ eV} = V \times e$  (volt x elementary charge (1))

- For example what is energy in eV in a circuit with a 20 volt electric field potential and

the charge flow of 40 electron

charges;  $E = 20V \times 40e = 800 \text{ eV}$

- $E_{(eV)} = V_{(V)} \times Q_{(C)} / 1.602176565 \times 10^{-19}$  (Volts to eV calculation with coulombs)
- To convert from eV to V with e an elementary charge of 1;  $V = eV / e$ : for example  $100 \text{ eV} / 1 = 100 \text{ V}$
- To convert from eV to V with coulombs,  $V_{(V)} = 1.602176565 \times 10^{-19} \times E_{(eV)} / Q_{(C)}$

#### Electron Charge e :

- charge carried by a single electron
- $e^- = \text{electron charge} = -1.602176565 \times 10^{-19} \text{ C/electron}$
- 1 coulomb (C) =  $6.242 \times 10^{18}$  electrons

#### Electron Gun:

Also known as an electron emitter source, a typically system produces a collimated electron beam that has a precise kinetic energy. A typical thermionic system is comprised of a heated cathode, electrodes to generate an electric field, including a Wehnelt electrode to initially focus the beam and an anode, with a large potential difference with the cathode to accelerate and further focus the electrons. The system typically needs to function in a moderate to high vacuum and requires several power supplies for various components of the systems.

#### Electron mass:

- $9.11 \times 10^{-31} \text{ kg/electron}$

#### Emission Current:

(1) For Electron Guns: The total electron current that leaves the cathode and goes to ground (in mA or  $\mu\text{A}$ ); equivalent to all the current that exits the gun (beam current) plus all the current that lands on grounded elements within the gun.

(2) For Ion Guns: The total ion current produced by the ion source that goes to ground (in mA or  $\mu\text{A}$ ), directly measured in only some gun; not identical to the electron emission current which causes gas ionization or to the ion beam current which exits from the gun; equivalent to all the current that exits the gun (beam current) plus all

---

the current that lands on grounded elements within the gun.

**Encoder Wheel:**

A disc-shaped rotary switch (thumb wheel) on early power supplies with FlexPanel control that adjusts the programming signal of a selected supply or menu item. The rate of increase/decrease varies with the speed of rotation.

**Emitter:**

A device for emitting charged particles, i.e. a cathode to emit electrons, or an anode to emit charged ions.

**Energy**

(1) Kinetic energy of a charged particle emitted from the gun; equal to  $\frac{1}{2}mv^2$ , where  $m$  is the particle mass and  $v$  is the particle velocity; this energy is achieved by applying accelerating voltages to elements of the gun. 1 eV is equivalent to the energy obtained by a particle with unit electronic charge ( $1.6021 \times 10^{-19}$  coulombs) accelerated by a 1 volt potential difference and is equal to  $1.6021 \times 10^{-19}$  joules.

(2) The energy power supply that provides the electron/ion accelerating voltage; other power supplies can be referenced to this power supply. Also, the potentiometer and switch that control the energy power supply.

(3) For Ion Guns: also called Ion Energy.

**Extract (also G-2):**

(1) The second electrostatic lens element in the ion gun, which provides the potential difference that extracts the ions from the ion source region and accelerates them; usually a tube with an aperture (hole); often grounded, but may be referenced to Energy and have a variable voltage.

(2) The power supply that applies voltage to the extract element in some ion guns. Also, the potentiometer that controls this power supply.

(3) For Ion Guns: also called Ion Energy

**Faraday cup:**

A particle-collecting device consisting of a shielded cup and output wire; usually used with an ammeter in series to ground to measure beam current.

**Feedthrough:**

Electrical, gas, mechanical and other connections (wires, tubing, insulators etc.), usually from vacuum to non-vacuum parts of a system.

**Filament, Fil+, Fil-:**

(1) The wire heating part of the standard cathode or ion source.

(2) Also the positive and negative leads to the cathode or ion source.

**Firing Unit:**

(1). For Electron Guns: The part of the gun that emits the electrons; the assembly that can be replaced as a unit, usually consisting of the complete triode (cathode, grid and anode) or just the cathode and grid; in some guns, only a part of the firing unit called the firing unit cartridge needs to be replaced.

(2) For Ion Guns: The part of the gun that produces the ions; the assembly that can be replaced as a unit, usually consisting of the gas ion source or alkali metal ion cartridge, the grid and sometimes extract elements.

**Flange Multiplexer:**

A vacuum component with one flange surface (usually 2.75" CF) to connect to a vacuum port and five smaller flanges (usually 1.33" CF) for various feedthrough and cable connections; used as the mounting system on many electron/ion guns.

**FlexPanel:**

(1) A Kimball Physics power supply control panel design featuring a digital screen,

---

selector buttons and a single variable control (up/down arrows).

(2) More specifically, the set of digital interface controller boards designed by Kimball Physics, used in power supplies for programmed control of supplies, metering, and display on a digital vacuum fluorescent screen. Power supplies with this controller are considered FlexPanel units even if they are “black box” style with no manual controls or display screen.

**Focus (also Lens):**

(1) One or more electrostatic or magnetic lens elements in the electron/ion gun that control the beam spread; usually controls spot size at the target; in some guns, may control beam divergence at points internal to the gun.

(2) The power supply that applies voltage or current to the focus/lens element(s). Also, the potentiometer that controls this power supply.

**Gaussian:**

A normal distribution curve: even if a beam current distribution is not strictly Gaussian, this term is used to describe distributions that are brightest in the center of the spot and fade off at the edges.

**G-1, G-2, G-3, G-4:**

The series of apertures and electrostatic lens elements in the electron/ion gun labeled by their position (first, second, third, fourth) instead of by common names.

**Grid (also G-1 or Wehnelt):**

The first electrostatic lens element in the electron/ion gun; the second element of the triode in electron guns; provides the first control and usually focuses the beam to a crossover; usually a tube with an aperture (hole), but may be a mesh (like a screen); can be used to cut-off the beam in most guns.

-The power supply applies voltage to the grid element. Also, the potentiometer that controls the grid power supply.

-This is all commonly referred to as the Wehnelt or grid cap.

**Gun Bias Voltage:**

This is the voltage between the filament and Wehnelt cap or grid that shapes the e- beam at its origin and keeps the filament current (feedback) at saturation.

**High Brightness:**

In an electron gun system, typically a cathode emitter that enables emission of electrons into a beam with a high current density.

**H.V.:**

High Voltage, sometimes used as a label on the Energy switch or to refer to an additional Power Supply unit needed for a high current electron gun.

**Insertion Length:**

The length of the gun that is in vacuum, from the vacuum sealing surface to the beam-exit end of the gun.

**Ion Cage:**

The part of the ion source where the gas is ionized in some ion guns, consisting of a wire coil that accelerates the ionizing electrons.

**Ion Emission Current:**

The total ion current produced by the ion source (in mA or  $\mu$ A); not identical to the electron emission current which causes ionization or to the ion beam current which exits from the gun.

**Ionization:**

(1). The production of a charged particle (ion) from a neutral atom or molecule by either the loss of one or more electrons (positive ion), or gain of one or more electrons (negative ion).

(2) Electron impact ionization- Ionization by bombarding a gas with electrons, which collide with and remove electrons from the neutral gas molecules producing positive ions.

(3) Surface ionization- Ionization by heating a solid alkali metal compound so that ions are produced in a solid-solid chemical reaction.

**Ion Source:**

---

(1) The structure in the ion gun that produces ions; for gas ions, usually consists of an electron-emitting filament, shields and an ion cage; for alkali metal ions, consists of a chemical cartridge and its heating filament.

(2) Another term for a complete ion gun; the entire unit made up of the ion source which creates the ions, and the electrostatic lens system that controls and focuses the ions to produce the final ion beam.

**Iridium:**

A conductive material used to mount or support materials used as electron emitters (Yttrium)

**Kelvin (K):**

- absolute temperature
- 0 Kelvin = absolute zero
- 0 Celsius (C) = 273.16 degrees (K) = triple point of water

**Lanthanum Hexaboride (LaB<sub>6</sub>):**

A cathode material with low work function and long lifetime; a single crystal structure cut and mounted on a heating rod; used as a cathode in some electron guns to produce electrons that can be focused to a small spot.

**LaB<sub>6</sub> Cathodes:**

Kimball Physics manufactures several cathodes based on LaB<sub>6</sub>, including single-crystal, guard ring, and planner disc.

- Kimball Physics references:

**LabVIEW™:**

National Instruments Inc. software used by Kimball Physics to develop programs to control the electron or ion gun power supply; allows the system to be operated remotely via a virtual instrument panel on a computer screen; also allows automated collection of data.

**Langmuir Equation:**

Used to determine brightness of thermionic cathode a is function of emitted current density, temperature, and beam energy.

**Lens (also Focus):**

(1) One or more electrostatic or magnetic lens elements in the electron/ion gun that control the beam spread; usually controls spot size at the target; in some guns, may control beam divergence at points internal to the gun.

(2) The power supply that applies voltage or current to the focus/lens element(s). Also, the potentiometer that controls this power supply.

**Perveance:**

A quantitative term used in the description of charged particle beams that indicates how significant the space charge effect is on the beams motion.

**Plasma:**

A state of matter in which outer electrons are not held to the rest of the atoms or molecules, so there are free electrons and free ions in a gas-like cloud; used as a source for ions in some guns.

**Power Supply:**

(1) The complete unit containing electronic circuitry, meters, and controls, that takes power input from the outside line and converts it to voltages and currents required to run the electron/ion gun.

(2) An individual group of electronic circuits, contained in the larger unit, that produces voltage/current applied to a specific element in the electron/ion gun.

**Pulsing (also Beam Pulsing):**

Rapidly stopping and starting the flow of electrons/ions, with a variable width (time on/off) and variable repetition rate; usually accomplished by rapidly switching the grid voltage to the cut-off potential; can be controlled in one of several ways: manual control with the



---

grid potentiometer, remote computer control, dual grid power supplies with a TTL switching signal, capacitive pulsing with a pulse junction box and an external pulse generator.

**Pulse Junction Box:**

A capacitor-containing device that combines input from an external pulse generator and the voltage from the grid power supply to vary the voltage applied to the grid; used to turn the beam current on and off rapidly; can be a separate box or part of a Grid Pulsing Cable.

**Rastering:**

Movement of the electron/ion beam in a synchronized pattern sweeping back and forth (X direction) while simultaneously moving down more slowly (Y direction), then returning to the start position without sweeping back and forth (retrace); used to cover an area of the target region fairly uniformly over time; produced by two synchronized cyclical voltages or currents applied to the X and Y deflection plates or deflection coils in the gun. Rastering may also be non-synchronized, in which case there is no retrace segment

**Refractory Metal Cathode:**

A cathode based on a class of metals that are extraordinarily resistant to heat and wear. Elements such as niobium, molybdenum, tantalum, tungsten and rhenium, with their melting points above 2000C, their high hardness and density, and relative chemical inertness are typically included in this group.

**Richardson Constant  $A_R$ :**

- $A_R = A/m^2K^2$
- $A_R$  (LaB<sub>6</sub>) = 29
- $A_R$  (CeB<sub>6</sub>) = 3.6

**Richardson-Dushman Equation:**

Equation used to estimate the current density (emission) from thermionic heated cathode electron source.

- $\sigma = j = A/m^2$
- $\sigma = j = A_R T^2 \exp(-e\Phi/kT)$
- $= A_R T^2 e(-e\Phi/kT)$
  
- $A_R$  = Richardson constant
- T = temperature K
- $\Phi$  = work function eV
- k = Boltzmann's Constant
- -e = electron charge =  $-1.60 \times 10^{-19}$   
C/electron

**Rocking (also Beam rocking):**

Movement of the electron/ion beam in an unsynchronized pattern sweeping back and forth (X direction), while simultaneously moving up and down more slowly (Y direction), with no retrace; used to cover an area of the target region uniformly over time; produced by magnetic fields from varying currents in electrical coils around the beam in the gun.

**Schottky Effect:**

The effect where a strong electric field (E) at the cathode surface enhances cathode emission current density and lowers the barrier to thermionic emission.

Secondary emission: a phenomenon whereby a primary incident particle of sufficient energy, hitting or passing through a surface, induces the emission of secondary particles.

**SEM:**

Scanning electron microscopy

**Source:**

(1) For Electron Guns: The power supply that produces the voltage and current that heats the

---

cathode causing it to emit electrons. Also the potentiometer and switch that control the source power supply.

(2) For Ion Guns: The power supply that produces the voltage and current that heats the filament of the ion source, either causing the filament to emit electrons for gas ionization, or heating the alkali metal cartridge for solid ionization, depending on gun. The potentiometer and switch that control the source power supply.

(3) The cable and power supply connector that include the source (filament), grid and other leads, depending on the electron/ion gun; the H.V. Multiconductor Source Cable.

(4) Ion source See Ion Source above.

**Space Charge:**

The electric field resulting from the presence of several charged particles (electrons and ions) in a region of space; can limit the emission of more electrons from a cathode or limit the ions extracted from an ion source; can also affect the path of charged particles through a region.

**Spot:**

The electron/ion beam at the target area; can be seen as a (usually circular) light area on a phosphor screen.

**Spot Size:**

The diameter (in mm or  $\mu\text{m}$ ) of a circle that contains most of the electron/ion beam at a given distance from the gun (working distance); measured either visually on a phosphor screen, or electrically by sampling beam current density across the spot with a Faraday cup; FWHM (full-width-half-max) measurement is the width that includes all beam current densities greater than half the maximum density.

**Sterad: Steradian:**

the SI unit of solid angle, equal to the angle at the center of a sphere subtended by a part of the surface equal in area to the square of the radius.

**Tantalum:**

A refractory metal commonly used as the electron emitting material in a cathode.

**Tantalum Disk Cathode:**

A Kimball Physics cathode electron emitter

- Kimball Physics resource:
- ES-042 Tantalum Disc Cathode
- ES-042 Tantalum Disc Cathode Tech Doc

**Target:**

The place (a plane) at which the electron/ion beam is aimed.

**TEM:**

Transmission electron microscope: a form of microscope where the image is derived from electrons passing through a sample, and where there is variable passage of electrons through the various spatial regions (x, y) of the specimen with the variable attenuation used to create an image. At lower magnifications TEM image contrast is due to differential absorption of electrons related to differences in material due in composition or thickness. For higher magnifications, more complex wave interactions occur and modulate the intensity of the image, requiring more advanced methods of images analysis.

**Tetrode:**

An electron extraction and acceleration unit used in some electron guns, consisting of four elements: cathode, grid, a variable first anode, and a grounded second anode.

**Thermionic Emission:**

Emission of electrons or ions from a metal or other coating material by heating the emitter in order to impart enough energy to overcome the work function of the emitting material.

**Triode:**

The basic electron extraction and acceleration unit of most electron guns, similar to a vacuum tube, consisting of three elements: cathode, grid (Wehnelt), and anode.

---

**Tungsten:**

A refractory metal commonly used in cathodes, because of its high melting point, either as the electron emission surface or as the heater leg.

**Vacuum:**

Typically measured in torr (760 torr = 1 atmosphere). Several types of pumping systems are used to lower pressure to create vacuum- roughing pumps, ion pumps and getter pumps.

- Low vacuum  $\sim 10^{-4}$  torr
- High vacuum  $\sim 10^{-7}$  torr
- Note that a high vacuum means lower pressure and less gas molecules.

**Vapor pressure:**

The pressure exerted by a vapor in equilibrium with its condensed phases (liquid or solid) in a closed space at a given temperature. Vapor pressure increases with temperature.

Equilibrium vapor pressure can be defined as the pressure reached when a condensed phase is in equilibrium with its own vapor. In the case of an equilibrium solid, such as a crystal, this can be defined as the pressure when the rate of sublimation (solid directly to gas) of a solid matches the rate of deposition of its vapor phase. For most solids this pressure is very low. However, with cathodes functioning at elevated temperature with small physical dimensions, the magnitude of physical structure loss secondary to sublimation is significant and limits the lifetime of the cathodes.

**Wehnelt:**

-Another term for grid, G1 or cap, from the physicist, A. Wehnelt.

-Creates a secondary electric field in the vicinity of the cathode to influence /control electron emission

**Work function:  $\Phi$ , W:**

A material-dependent measure of the work per unit charge needed to emit a charged particle (usually an electron) from the emitting material into the vacuum just adjacent to the material surface.

- $\Phi$  (LaB<sub>6</sub>) = 2.70 eV
- $\Phi$  (CeB<sub>6</sub>) = 2.65 eV
- $\Phi$  (Tungsten) = 4.5 eV

**X axis, Y axis:**

Arbitrarily-defined horizontal and vertical lines used as reference for motion of the beam in the target area for deflection, rastering, or alignment; correspond to the position of four deflection plates in the gun.

**Yttrium:**

Material with atomic number of 39 that is a transition metal used for fabrication of electron sources. At Kimball Physics, the oxide of Yttrium, Y<sub>2</sub>O<sub>3</sub>, is used as a coating on the surface of Iridium and used as an electron emission source.

**Yttria-Coated Iridium Cathodes:**

A Kimball Physics thermionic cathode that is designed for more rugged applications and is more resistant to oxidation in less favorable vacuum conditions.

**Z axis:**

A reference line through the exact center of the electron or ion gun, in the direction of electron or ion travel; the approximate beam line.

**Zoom Lens:**

A series of electrostatic lens elements in the electron/ion gun (including anode or G-2, focus, and sometimes other lenses) that both focuses and accelerates / decelerates the beam; allows spot size and beam current to remain roughly constant as energy is varied.

---

Notes:

1. Charts /graphs show typical performance, data is for guidance only
2. It is not necessarily possible to achieve all maximum specifications simultaneously.
3. Specifications Subject to Change Without Notice.
4. DE Altobelli 6/6/2023
5. Beta Version New Document- Recommendations, revisions, corrections are welcomed.

Document: Electron\_Gun\_Tutorial\_Resource\_2023\_0606  
COPYRIGHT KIMBALL PHYSICS 2023, ALL RIGHTS RESERVED